

Gordon Kearney¹, Thorsten Bernsmann²
¹Waters Corporation, Manchester, UK. ²CVUA Munster, 48151 Munster, Germany

INTRODUCTION

Under poor storage conditions, agricultural produce is often subject to fungal contamination. *Penicillium* and *Aspergillus* molds, which produce the natural product ochratoxin A, may be found on stored coffee beans, grapes and grain¹. Ochratoxin A is carcinogenic in rats and may be associated with human kidney disease. Because of this a Maximum Residue Level (MRL) in grain of 5 µg/Kg has been proposed in the European Union². The structure of ochratoxin A is shown in Figure 1.

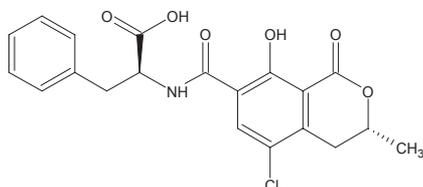


Figure 1. Ochratoxin A.

Many other foodstuffs, including beer and wine, may also contain unacceptably high levels of ochratoxin A. For this reason, a rapid and simple method was developed for its determination in red wine. No extraction or cleanup was performed. An aliquot of wine was analysed directly and a fast chromatographic gradient allowed a sample to be analysed every 10 minutes.

METHOD

All experiments were conducted using a Waters® Alliance® HPLC system, an XTerra® 2.1 mm by 30 mm HPLC column and a Waters Micromass® Quattro micro™ API triple quadrupole mass spectrometer. Mobile phase A was water with 1% formic acid. Mobile phase B was acetonitrile with 1% formic acid. 50 µL of red wine was loaded onto the HPLC column and the following solvent gradient was run at a flow rate of 0.4 mL/min.

Time	0 mins	5% B
Time	1 min	5% B
Time	3 min	100% B
Time	6 min	100% B

The eluent was interfaced to the electrospray source of the mass spectrometer in positive ion mode. The MS/MS transitions selected for quantification and confirmation of the compound are shown in Table 1.

Precursor m/z	Product m/z	Dwell Time (s)	Cone Voltage (V)	Collision Energy (eV)
403.9	238.9	0.1	21	25
403.9	357.9	0.1	21	15
405.9	240.9	0.1	21	25

Table 1. MS/MS conditions for ochratoxin A.

In order to avoid contamination of the ion source, the LC eluent was diverted to waste for the first three minutes of each run.



Waters Micromass Quattro micro API triple quadrupole mass spectrometer.

RESULTS

Figure 2 shows the full scan MS spectrum of ochratoxin A under positive electrospray conditions.

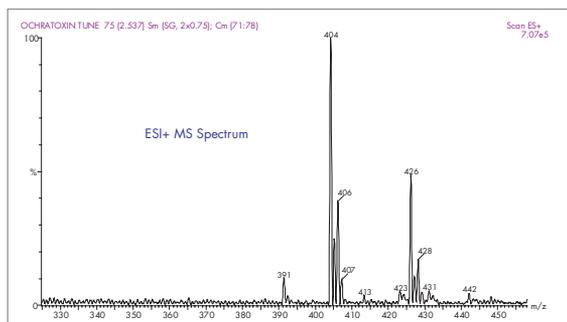


Figure 2. MS spectrum of ochratoxin A.

Figure 3 shows the chromatograms obtained from the analysis of red wine spiked at 0.25 µg/L.

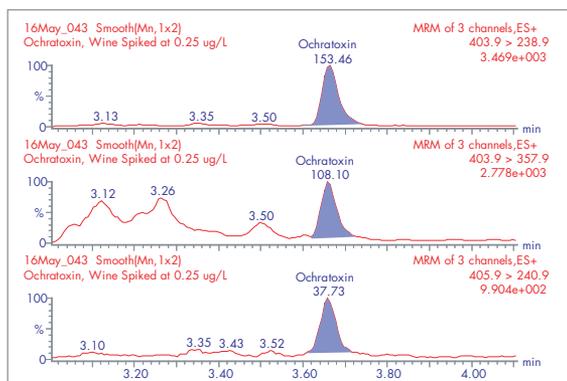


Figure 3. Chromatograms for 0.25 µg/L ochratoxin A in red wine.

A calibration graph was generated between 0.024 and 2.39 µg/L and is shown in Figure 4.

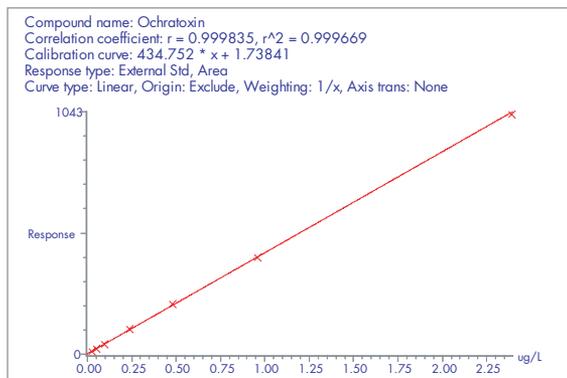


Figure 4. Calibration graph for ochratoxin A.

CONCLUSION

Given the concerns over the potential toxic effects of ochratoxin A in the human diet, it is necessary to study its occurrence in various foods. The method described above allows for the rapid and sensitive analysis of this compound in red wine, without the need for extraction and cleanup procedures.

REFERENCES

1. van Egmond HP, Speijers GJA. J. Nat. Toxins, 1994, 3: 125.
2. Krivobok S, Seigle-Murandi F, Steiman R, Creppy EE. System. Appl. Microbiol. 1995; 18: 455.

Sales Offices

AUSTRIA AND EXPORT (CENTRAL EUROPE, CIS, MIDDLE EAST,
INDIA AND INDIA SUBCONTINENT)
43 1 877 18 07

AUSTRALIA 61 2 9933 1777

BELGIUM AND LUXEMBOURG
32 2 726 1000

BRAZIL 55 11 5543 7788

CANADA 800 252 4752 X2205

CZECH REPUBLIC 420 2 617 11384

DENMARK 45 46 59 8080

FINLAND 358 09 506 4140

FRANCE 33 1 3048 7200

GERMANY 49 6196 400600

HONG KONG 852 29 64 1800

HUNGARY 36 1 350 5086

INDIA 91 80 837 1900

IRELAND 353 1 448 1500

ITALY 39 02 27 4211

JAPAN 81 3 3471 7191

KOREA 82 2 3284 1300

MEXICO 52 55 5524 7636

THE NETHERLANDS 31 76 508 7200

NORWAY 47 6 384 6050

PEOPLE'S REPUBLIC OF CHINA
86 10 8451 8918

POLAND 48 22 833 4400

PUERTO RICO 787 747 8445

RUSSIA/CIS 7 095 931 9193

SINGAPORE 65 6278 7997

SPAIN 34 93 600 9300

SWEDEN 46 8 555 11 500

SWITZERLAND 41 62 889 2030

TAIWAN 886 2 2543 1898

UK 44 208 238 6100

U.S.A. AND ALL OTHER COUNTRIES:

WATERS CORPORATION
34 Maple St.
Milford, MA 01757 U.S.A.
T: 508 478 2000
F: 508 872 1990

WATERS CORPORATION

34 Maple St.

Milford, MA 01757 U.S.A.

T: 508 478 2000

F: 508 872 1990

www.waters.com

Waters
RIGHT ON TIME.



Waters, Micromass, XTerra, Quatro micro and Alliance are trademarks of Waters Corporation.
All other trademarks are the property of their respective owners.
©2003 Waters Corporation December 2003 720000792EN DE-PDF
Produced in the United Kingdom.

