

Ashley Sage
¹Waters Corporation, Manchester UK

INTRODUCTION

Why Exact Mass?

As with every analytical tool, the more accurate your measurement the more confident you can be with your results. For example, when carrying out a metabolism investigation following an *in vivo* experiment are you sure that a peak detected at m/z 342.1 is an actual metabolite or a matrix related component or even a solvent contaminant? Mass measuring the peak of interest to 4 decimal places will give you that confidence.

By reducing the error on a mass measurement, the possible elemental formulae that can fit a particular measurement significantly reduce, therefore providing better specificity. Figure 1 clearly shows this. Such mass measurement accuracy is not routinely available with typical scanning instruments such as quadrupole or ion trap mass spectrometers. For routine and easy-to-use exact mass measurement, orthogonal acceleration time-of-flight (oa-TOF) mass spectrometry provides the solution.

Exact mass measurement on an oa-TOF can help to:

- Determine the elemental compositions of analytes.
- Distinguish between nominally isobaric components.
- Distinguish between analyte and matrix related material.
- Improve database search specificity.
- Increase your confidence in analyte identification.

For LC/MS applications, the Waters® Micromass® LCT Premier™ benchtop oa-TOF mass spectrometer provides the user with the ability to acquire high resolution MS data and thus generate an exact mass measurement. In addition to the next generation of instrument electronics and oa-TOF optics, LCT Premier is equipped with LockSpray™ as standard. LockSpray is a novel dual electrospray ion source that was designed for dedicated exact mass measurement. Your LC column is connected to the analyte sprayer while the single point reference mass, used to exact mass measure your data, is introduced into the reference sprayer.

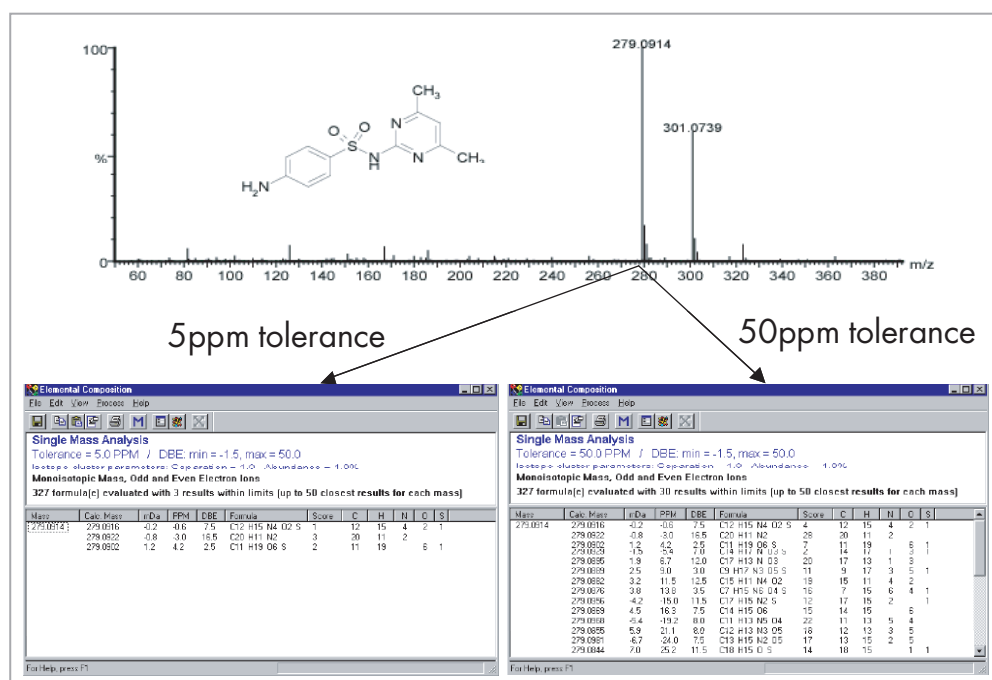


Figure 1. Improving Specificity Through Exact Mass Measurement.

The eluent from each of the two sprayers is kept completely separate by a baffle that is indexed via software control for total data integrity. During an experiment, data from the reference sprayer is typically sampled once every 10 secs and the reference data is automatically applied to the analyte data to provide an exact mass measurement. LockSpray overcomes issues associated with traditional exact mass measurement, i.e. post-column addition of a lock mass and potential ionization interferences between analytes and lock mass etc.

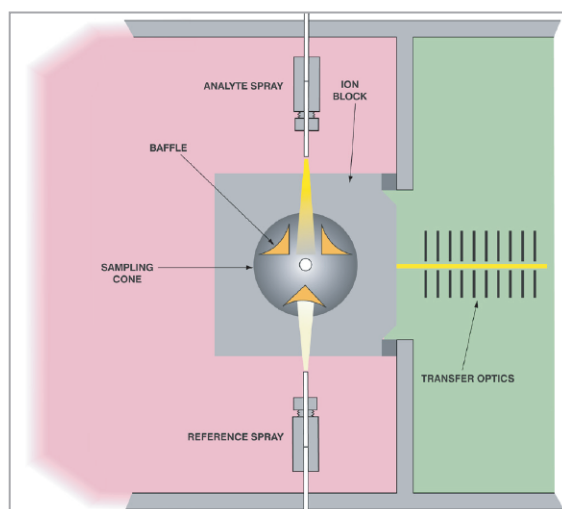


Figure 2. LockSpray in Analyte Position.

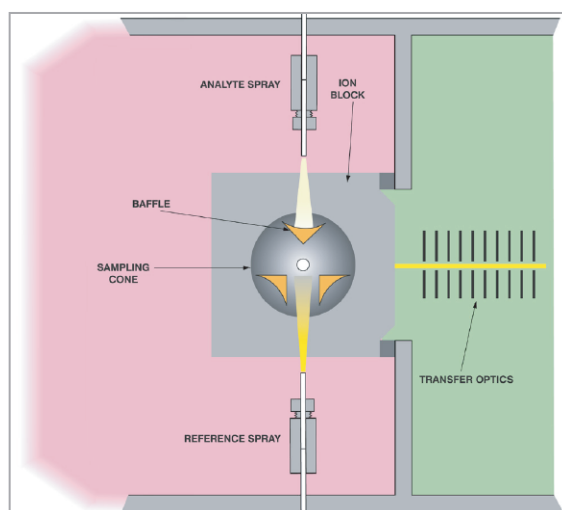


Figure 3. LockSpray in Reference Position.

Examples of Mass Measurement Accuracy on LCT Premier

1) Single Analyte Exact Mass Measurement.

- LockSpray with positive ion electrospray was used.
- Raffinose (500 pg/ μL in 50/50 MeCN/ H_2O) was infused into the analyte sprayer over 30 mins at 5 $\mu\text{L}/\text{min}$. The $[\text{M}+\text{Na}]^+$ ion was measured ($\text{C}_{18}\text{H}_{31}\text{O}_{16}\text{Na}$, m/z 527.1588).
- Leucine enkephalin (50 pg/ μL in 50/50 MeCN/ H_2O + 0.1% formic acid) was infused into the reference sprayer at 5 $\mu\text{L}/\text{min}$. The $[\text{M}+\text{H}]^+$ ion was used as the lock mass ($\text{C}_{28}\text{H}_{37}\text{N}_5\text{O}_7$, m/z 556.2771).
- The reference sprayer was sampled every 10 secs and 10 measurements were taken over the 30 min acquisition.
- The spectral resolution was 10,000 FWHM.

Table 1 below lists the exact mass measurements along with the calculated mDa and ppm errors from the theoretical. The RMS ppm error for all 10 measurements was calculated at 1.32 ppm.

| Actual Mass | Measured Mass | mDa Error | ppm Error |
|---------------|---------------|-----------|-----------|
| 527.1588 | 527.158 | -0.8 | -1.52 |
| 527.1588 | 527.1581 | -0.7 | -1.33 |
| 527.1588 | 527.1588 | 0 | 0.00 |
| 527.1588 | 527.1578 | -1 | -1.90 |
| 527.1588 | 527.1581 | -0.7 | -1.33 |
| 527.1588 | 527.1578 | -1 | -1.90 |
| 527.1588 | 527.1582 | -0.6 | -1.14 |
| 527.1588 | 527.1595 | 0.7 | 1.33 |
| 527.1588 | 527.1583 | -0.5 | -0.95 |
| 527.1588 | 527.1585 | -0.3 | -0.57 |
| RMS ppm error | | | 1.32 |

Table 1. Exact Mass Measurement Results for Raffinose.

2) Multiple Analyte Exact Mass Measurement during an LC/MS Experiment.

- LockSpray with positive ion electrospray was used.
- An 8 component analyte mixture was separated using the following conditions:
 - HPLC: Waters Alliance® HT 2795.
 - Mobile phase A : H₂O + 0.1% formic acid.
 - Mobile phase B : MeCN + 0.1% formic acid.
 - Gradient : 95% A to 5% A over 10 mins.
 - Column : Waters Atlantis™ dC18, 2.1 x 100 mm, 3 μ .
 - Flow rate : 0.3 mL/min.
 - Injection volume : 10 μ L.
- Leucine enkephalin (50 pg/ μ L in 50/50 MeCN/H₂O + 0.1% formic acid) was infused into the reference sprayer at 5 μ L/min. The [M+H]⁺ ion was used as the lock mass (C₂₈H₃₇N₅O₇, m/z 556.2771).
- The reference sprayer was sampled every 10 secs.

Figure 4 shows a typical chromatogram obtained with the retention time for all 8 components. Table 2 lists the exact mass measurements obtained along with the calculated mDa and ppm errors from the theoretical. The RMS ppm error for all measurements was calculated at 1.8 ppm.

This example shows that the LCT Premier provides excellent mass measurement accuracy across a wide range of compounds with varying mass, physical properties and retention time during a reversed phase LC gradient. This provides improved specificity during an LC/MS experiment providing the analyst with increased confidence in the analytical result.

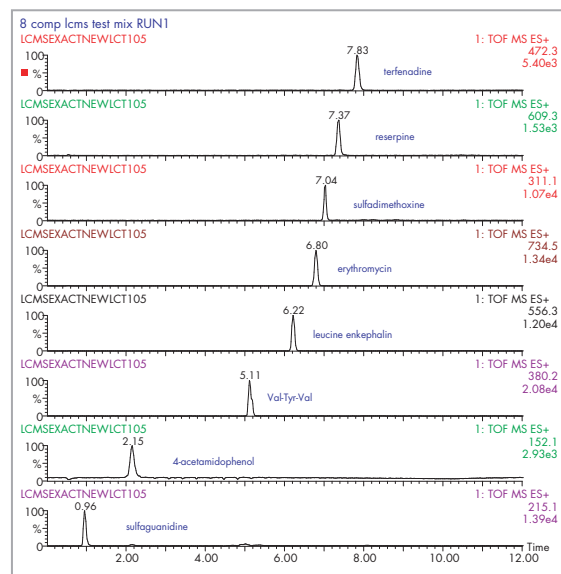


Figure 4. Retention Times for all 8 Components.

| Compound | Actual Mass | Measured Mass | mDa Error | ppm Error |
|--------------------|-------------|---------------|-----------|-----------|
| 4-acetamidophenol | 152.0711 | 152.0707 | -0.4 | -2.6 |
| sulfaguanidine | 215.0602 | 215.0597 | -0.5 | -2.3 |
| Sulfadimethoxine | 311.0814 | 311.081 | -0.4 | -1.3 |
| Val-Tyr-Val | 380.2185 | 380.219 | 0.5 | 1.3 |
| Terfenadine | 472.3215 | 472.321 | -0.5 | -1.1 |
| Leucine enkephalin | 556.2771 | 556.2775 | 0.4 | 0.7 |
| Reserpine | 609.2812 | 609.2828 | 1.6 | 2.6 |
| Erythromycin | 734.469 | 734.4695 | 0.5 | 0.7 |
| RMS ppm error | | | | 1.8 |

Table 2. Exact Mass Measurements of 8 Component Mix.

3) Exact Mass Measurement during Positive/Negative Ion Mode Switching.

Having the ability to carry out ionization mode switching in real time provides the analyst with an additional piece of information during the same LC/MS experiment. Ionization mode switching is often regarded as being useful for when the expected mode of ionization for a particular compound is unknown e.g. compound screening experiments. With the

additional function of being able to generate an exact mass measurement during ion mode switching experiment in real time, further confidence in that the correct ion has been detected is obtained.

With significant improvements in electronics and detector hardware, the Waters Micromass LCT Premier now has the possibility to switch ionization mode during a real time LC/MS experiment in 300 msec—providing an exact mass measurement and a sufficient number data points to define a typical chromatographic peak. The previous oa-TOF instrument, the LCT, was only able to switch ionization mode during a real time experiment in 1 second, due to the stabilization time required when the older generation high voltage electronics were switched polarity.

- LockSpray with positive/negative ionization mode switching was used.
- An Inter Scan Delay (ISD) of 0.3 secs was used.
- The reference sprayer was sampled every 5 secs.
- Raffinose ($C_{18}H_{32}O_{16}$, 500 pg/ μ L in 50/50 MeCN/ H_2O) was infused into the analyte sprayer at 5 μ L/min.
 - $[M+Na]^+ = 527.1588$ m/z.
 - $[M-H]^- = 503.1612$ m/z.
- Leucine enkephalin ($C_{28}H_{37}N_5O_7$, 50 pg/ μ L in 50/50 MeCN/ H_2O + 0.1% formic acid) was infused into the reference sprayer at 5 μ L/min.
 - $[M+H]^+ = 556.2771$ m/z.
 - $[M-H]^- = 554.2615$ m/z.

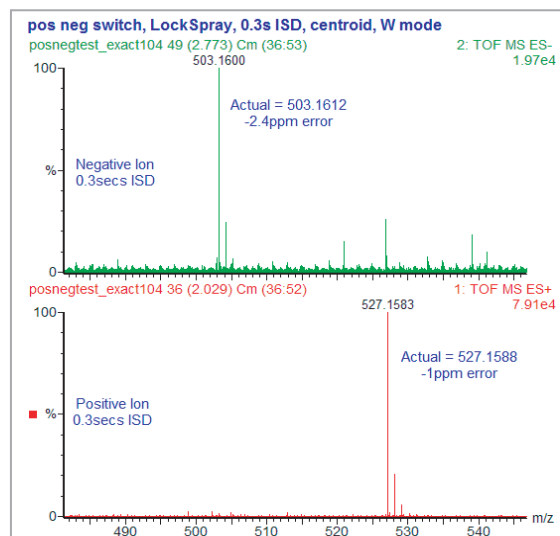


Figure 5. Positive/Negative Ion Mode Switching with LockSpray.

Figure 5 shows a typical positive and negative ion spectrum for raffinose that was acquired during the experiment. The exact mass measurements shown are representative of the data that is typically obtained during a positive/negative ion mode switching experiment with LCT Premier.

CONCLUSION

For LC/MS applications, the LCT Premier provides the user with the ability to acquire high resolution MS data and therefore the ability to generate an exact mass measurement to within 3 ppm. Using the LockSpray ion source, traditional problems associated with exact mass measurement during LC/MS acquisitions are overcome—no multi point reference masses, no post-column addition of the lock mass, no ionization interference between your analytes and lockmass, and no nominal mass interferences. Having the ability to acquire an exact mass measurement aids determination of elemental composition for a particular analyte of interest, which provides the analyst a high degree of specificity and therefore a high level of confidence in the analytical result.

LCT Premier—exact mass measurement made easy for LC/MS applications.

Sales Offices

AUSTRIA 43 1 877 18 07

AUSTRALIA 61 2 9933 1777

BELGIUM AND LUXEMBOURG
32 2 726 1000

BRAZIL 55 11 5543 7788

CANADA 800 252 4752 X2205

CZECH REPUBLIC 420 2 617 11384

DENMARK 45 46 59 8080

FINLAND 358 9 506 4140

FRANCE 33 1 3048 7200

GERMANY 49 6196 400600

HONG KONG 852 29 64 1800

HUNGARY 36 1 350 5086

INDIA 91 80 837 1900

IRELAND 353 1 448 1500

ITALY 39 02 27 4211

JAPAN 81 3 3471 7191

KOREA 82 2 3284 1300

MEXICO 52 55 5524 7636

THE NETHERLANDS 31 76 508 7200

NORWAY 47 6 384 6050

PEOPLE'S REPUBLIC OF CHINA
86 10 8451 8918

POLAND 48 22 833 4400

PUERTO RICO 787 747 8445

RUSSIA/CIS 7 095 931 9193

SINGAPORE 65 6278 7997

SPAIN 34 93 600 9300

SWEDEN 46 8 555 11 500

SWITZERLAND 41 62 889 2030

TAIWAN 886 2 2543 1898

UK 44 208 238 6100

U.S.A. AND ALL OTHER COUNTRIES:

WATERS CORPORATION

34 Maple St.

Milford, MA 01757 U.S.A.

T: 508 478 2000

F: 508 872 1990

www.waters.com

WATERS CORPORATION

34 Maple St.

Milford, MA 01757 U.S.A.

T: 508 478 2000

F: 508 872 1990

www.waters.com

Waters

Opt-In! My Profile
www.waters.com

Waters, Micromass, LCT Premier, LockSpray, Alliance and Atlantis are trademarks of Waters Corporation.

All other trademarks are the property of their respective owners.

©2004 Waters Corporation April 2004 720000810EN DE&LW-UL

