

Mike McCullagh and Tabisam Khan, Waters Corporation, Manchester, UK

Objective

The aim of the study presented is to determine the presence of target Ginkgolides in Ginkgo biloba leaf extract. The data presented shows Ginkgolides determined to be present in the extract analyzed. Using negative ion mode electrospray very good mass accuracy is achieved routinely with real time centroid data acquisition and lockmass correction. Acquisition of UV data in parallel is performed. The system used was comprised of a Waters® Alliance® HT 2795 Separations Module, 2996 PDA Detector, Symmetry® C₁₈ Column and Waters Micromass® LCT Premier™ oa-TOF Mass Spectrometer.

Introduction

Medical treatment has long been connected with the natural products and actives extracted from plants. This has led to many advances in medicines and has long been the pivotal force in fighting ailments. It has been known and appreciated for many years that certain plants can exhibit healing effects when prepared as part of a mixture, but when taken in isolation can cause detrimental effects. This knowledge is now being used more and more in sophisticated biological and bioanalytical studies of natural products. Despite advances both in utilizing synthetic approaches to drug design, and in sophisticated structure-activity studies, there is still a great need for compounds with a unique mechanism of action. Major breakthroughs have resulted primarily from the study of natural products. Some of the most important drugs have been isolated from plant sources; for example most antibiotics and anti-cancer drugs. The maidenhair tree, Ginkgo biloba, is an ancient Chinese plant that has been cultivated for its health promoting properties. Ginkgolides, the main active ingredients of Ginkgo biloba, have not only helped to explain the pharmacological basis of several traditional medicines, but have also

provided a valuable new class of therapeutic agents. Research on the biochemical effects of Ginkgo biloba extracts is still at a very early stage. Although the terpene fraction of Ginkgo biloba, which contains the ginkgolides, may contribute to the neuroprotective properties of the Ginkgo biloba leaf, it is also likely that the flavonoid fraction, containing free radical scavengers, is important in this respect. The structures of the ginkgolides are shown in Figure 4.

The LCT Premier is a newly developed benchtop oa-TOF (orthogonal acceleration time of flight) mass spectrometer. New hardware and software control technology has been incorporated to meet the increased analytical demands in the pharmaceutical, environmental and clinical applications arenas. The high duty cycle of TOF is utilized for qualitative studies, generating full spectra at high mass accuracy (<3 ppm RMS). The highly specific data generated provides an extra degree of information that aids interpretation of the data. Using real time exact mass centroid data acquisition with the negative ion electrospray, the extract of Ginkgo biloba leaf has been analyzed using the new LCT Premier oa-TOF. The LCT Premier oa-TOF is presented in Figure 1 and the schematic of the LCT Premier with the analyzer in W geometry is illustrated in Figure 2.



Figure 1. LCT Premier incorporating integral LockSpray™.

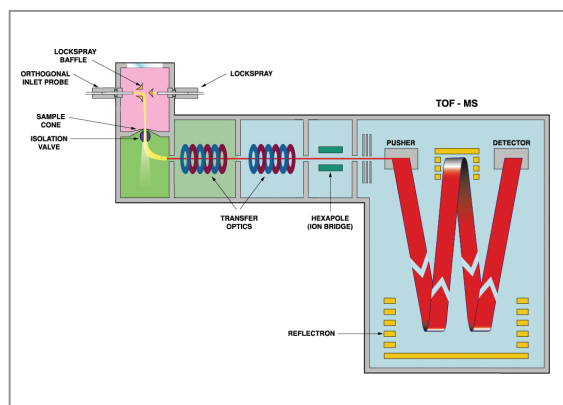


Figure 2. Oa-TOF schematic (W mode > 10000 FWHM).

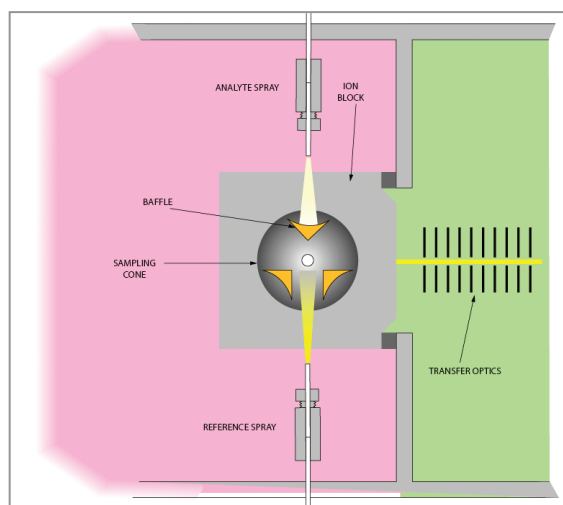


Figure 3. Schematic of the LCT Premier LockSpray source.

Experimental

Mass Spectrometer: Micromass LCT Premier oa-TOF

HPLC: Waters Alliance HT 2795 Separations Module

HPLC

Column: Waters Symmetry C₁₈ (250 mm x 4.6 mm, 5 µm particles) with guard column (20 mm x 3.9 mm, 5 µm particles)

Column temperature: 35 °C

Flow: 1 mL/min - split 1:4.

Mobile phase: A: H₂O (0.2% HCOOH), B: MeCN

Gradient: 0-10 min: 15% B; 10-40 min: 15-30% B; 40-50 min: 30-15% B

MS

Ionization mode: ESI Voltage -ve = 2.5kV.

Sample cone voltage: 100V.

Reference mass: Leucine enkephalin,
[M-H]⁻ = 554.2615

Acquisition parameters:

100 - 800 Da	1 spectrum/second
10500 FWHM	0.1 second inter scan delay

Extraction Procedure

The extraction procedure used by the China Pharmaceutical Company is described as follows;

- 50g Ginkgo Biloba leaf powder was refluxed with 200 mL 60% acetone for 1hr, then filtered.
- Extraction solvent was recovered and remaining residue was extracted with ethyl acetate.
- Ethyl acetate was recovered and remaining residue was extracted with water.
- Water was extracted with diethyl ether.
- Acetone, ethyl acetate and diethyl ether fractions were combined and evaporated.
- Samples were reconstituted in 60% acetone.

RESULTS

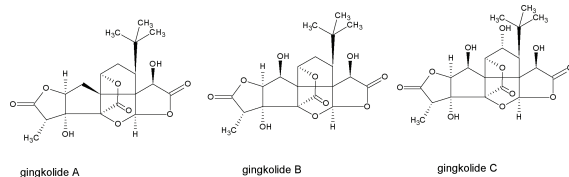


Figure 4. Chemical Structure of Ginkgolides.

The extract of *Ginkgo biloba* was analyzed using the LCT Premier in negative ion mode. Presented in Figure 4 are the structures of the Ginkgolides of interest. The negative mode total ion chromatogram obtained from the extract of Nanjing *Ginkgo Biloba* leaf is shown in Figure 5, where Ginkgolides A, B and C determined to be present are shown. From the chromatogram it can be seen that five major components were present in the purified extract. Under the chromatographic conditions used it was not possible to achieve complete separation of the Ginkgolides.

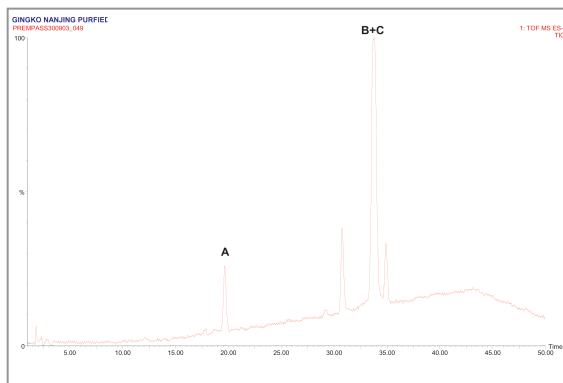


Figure 5. Negative mode total ion chromatogram obtained for the extract of Nanjing Ginkgolide *Biloba*.

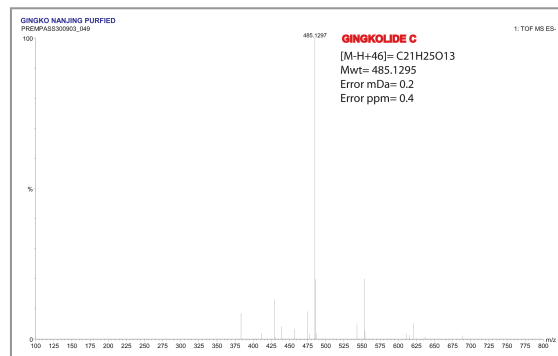


Figure 6. Exact mass spectrum obtained for Ginkgolide A.

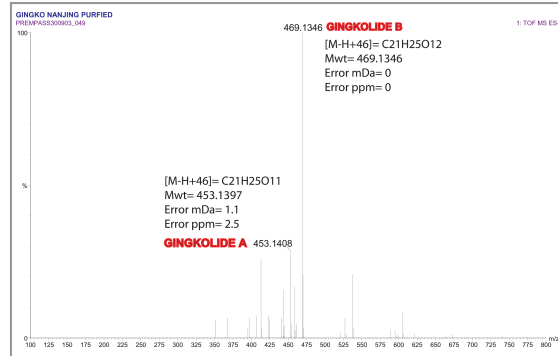


Figure 7. Exact mass spectrum obtained for Ginkgolide B and C.

The reference compound Leucine enkephalin ionizes in negative mode to produce $[M-H]^- = 554.2615$, and is sampled independently from the analyte spray to provide a lockmass. As shown in Figures 6 and 7 it is not seen in the analyte mass spectrum. Lockmass correction takes place automatically in real time, the independent sampling enhances the mass accuracy obtained. The exact mass spectrum obtained for Ginkgolide C is presented in Figure 6. Ginkgolide C was mass measured within less than 1 ppm error. For Ginkgolides A and B the exact mass spectrum is shown in Figure 7, where mass measurement has been obtained within 3 ppm error. For the Ginkgolides identified, in each case a formate adduct has been observed.

Conclusion

- Extracts of Gingko Biloba samples have been successfully analyzed allowing the specific Ginkgolide profile to be obtained.
- Analysis has been performed using centroid acquisition mode and exact mass measurement performed in real time.
- 0.4 ppm error was obtained for negative mode acquisition of Ginkgolide A.
- 0.0 ppm error was obtained for negative mode acquisition of Ginkgolide B.
- 2.5-ppm error was obtained for negative mode acquisition of Ginkgolide C.
- The LCT Premier oa-TOF with integral LockSpray and independent reference mass acquisition enables the routine rapid acquisition of highly specific data.
- Analyte confirmation is achieved using exact mass measurement and enabled complete confidence in the identification of the Ginkgolides.

Sales Offices:**AUSTRIA** 43 1 877 18 07**AUSTRALIA** 61 2 9933 1777**BELGIUM AND LUXEMBOURG**
32 2 726 1000**BRAZIL** 55 11 5543 7788**CANADA** 800 252 4752 X2205**CZECH REPUBLIC** 420 2 617 11384**DENMARK** 45 46 59 8080**FINLAND** 358 9 506 4140**FRANCE** 33 1 3048 7200**GERMANY** 49 6196 400600**HONG KONG** 852 29 64 1800**HUNGARY** 36 1 350 5086**INDIA** 91 80 2837 1900**IRELAND** 353 1 448 1500**ITALY** 39 02 27 4211**JAPAN** 81 3 3471 7191**KOREA** 82 2 820 2700**MEXICO** 52 55 5524 7636**THE NETHERLANDS** 31 76 508 7200**NORWAY** 47 6 384 6050**PEOPLE'S REPUBLIC OF CHINA**
86 10 8451 8918**POLAND** 48 22 833 4400**PUERTO RICO** 787 747 8445**RUSSIA/CIS** 7 095 931 9193**SINGAPORE** 65 6278 7997**SPAIN** 34 93 600 9300**SWEDEN** 46 8 555 11 500**SWITZERLAND** 41 62 889 2030**TAIWAN** 886 2 2543 1898**UK** 44 208 238 6100**US** 800 252 4752

Waters

WATERS CORPORATION

34 Maple St.

Milford, MA 01757 U.S.A.

T: 508 478 2000

F: 508 872 1990

www.waters.com

For Complete Confidence

Waters, Alliance, Symmetry, LCT Premier, LockSpray and Openlynx are trademarks of Waters Corporation.

All other trademarks are the property of their respective owners.

©2004 Waters Corporation Produced in the U.S.A. November 2004 720000980EN MC-PDF



001