

THE APPLICATION OF GC/MS FOR THE METABONOMIC ANALYSIS OF RAT URINE

Jennifer H. Granger¹, Doug Stevens², Nicholas Ellor², Robert S. Plumb¹, and Ian D. Wilson³

¹Waters Corporation, Milford, MA, USA , ²Waters Corporation, Beverly, MA, USA

³AstraZeneca, Alderley Park, Macclesfield, UK

INTRODUCTION

Metabonomics is the study of changes in the endogenous metabolites in biological fluids such as urine and plasma as a result of a toxic insult, disease state, or genetic difference. In order to detect these changes, information-rich techniques such as mass spectrometry coupled with a liquid or gas chromatographic separation process are employed. The greater the sensitivity and resolving power of the system, the more analytes detected and hence a greater chance of finding a biomarker of disease or toxicity. Capillary GC coupled to time-of-flight mass spectrometry, while requiring sample pre-treatment via derivatization, is relatively simple to use and provides the highly sensitive analyses required for multivariate statistical processing, as with Principal Components Analysis (PCA). Here we show how GC/MS(TOF) using the Waters Micromass® GCT™ Mass Detector has been employed with both EI and CI detection for the analysis of rat urine in a metabonomics study.

EXPERIMENTAL CONDITIONS

GC Conditions

Column: DB5MS 30 m x .25 mm ID x .25 µm film
Flow Rate: 1.5 mL/min. Helium
Injection Volume: 1 µL, split 10:1
Injector Temp: 275 °C
Temp Program: 50 °C for 4 min., ramp to 300 °C at 15 °C/min., hold for 5 min.
Solvent Delay: 6 min.

MS Conditions

Acquisition Range: 50 - 700 Da EI, 100 - 700 Da CI
Scan Time: 0.15 s
Interscan Delay: 0.05 s
Ionization: EI⁺, CI⁺ using ammonia
Lock Mass: tris-trifluoromethyltriazine (metri)
Source Temp: 180 °C EI, 175 °C CI
Transfer Line Temp: 250 °C

Animal Samples

Urine samples were collected from male and female obese (fa/fa) Zucker rats (BABU, Alderley Park; n = 10 of each sex; males age 13 weeks and females age 8 weeks) at two time periods, morning (AM) and evening (PM) to give a total of 40 samples. Samples were stored at -20 °C prior to analysis..

Sample Preparation

The urine samples were thawed and a 50 µL aliquot of each sample evaporated to dryness. A 250 µL aliquot of pyridine was added to each. This was immediately followed by the addition of 50 µL of BSTFA with 1% TCMS. Samples were then vortexed for 15 seconds and then heated to 80 °C for one hour. Each 300 µL derivatized sample aliquot was transferred to an autosampler vial with a 400 µL low-volume insert. Each sample was analyzed in triplicate.

RESULTS AND DISCUSSION

It is projected that type II diabetes will be the biggest killer in the western world by the year 2020. Thus, the biomarker information that can be derived from the study of Zucker rats (a naturally diabetic genetic strain) may hold important clues to controlling the disease.

Successful biomarker discovery is dependent upon high quality analytical data. The capillary gas chromatographic analysis of the rat urine samples produces a very information-rich complex EI chromatogram, generated by the GCT. The GCT is a time-of-flight (TOF) mass detector optimized for high resolution (7000 FWHM) and sensitivity for exact mass measurement. When operated in EI mode, the system is capable of generating library-searchable spectra with a high mass accuracy (<5 ppm error). In the event that an analyte is not contained in a library search, the elemental composition deduction capabilities of the GCT assist in the successful determination of unknowns.

Application NOTE

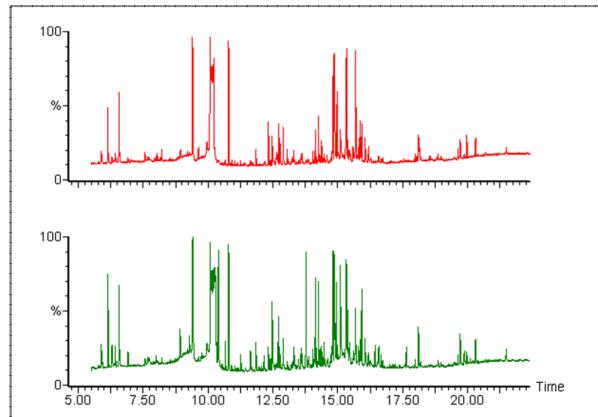


Figure 1.

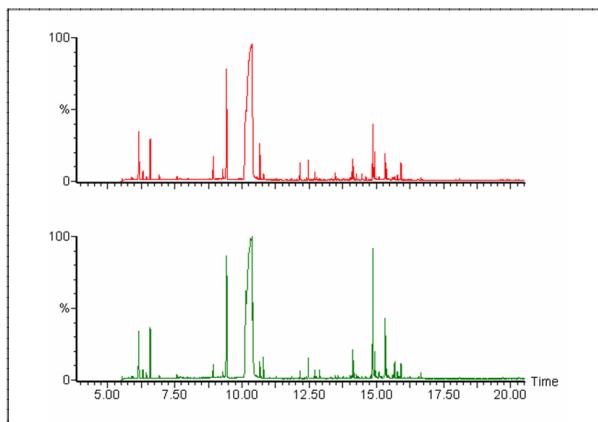


Figure 2.

A representative chromatogram from each of the Zucker and control animal groups is shown in Figure 1. While there are some subtle qualitative differences (e.g., the intensity of the peak at retention time 8.75 min.), the chromatograms are very similar. The CI GC/MS chromatogram from the same two samples is given in Figure 2. As we can see from the two data sets, the EI chromatogram is more information-rich than the CI analysis; however, the CI data provides molecular weight information which is not present in the EI data. The average peak width at the base for the EI GC/MS data was 0.05 min. with a peak capacity of approximately 400 for the twenty-minute separation.



Figure 3.

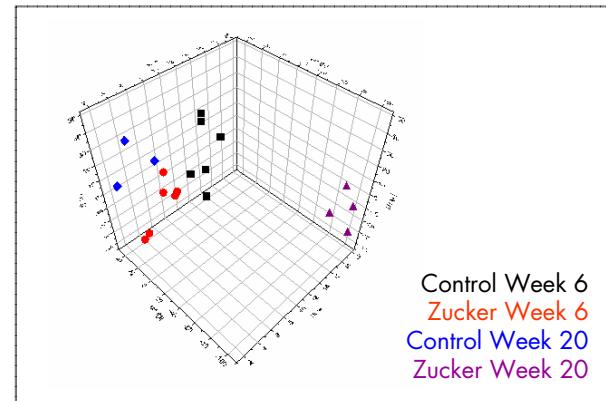


Figure 4.

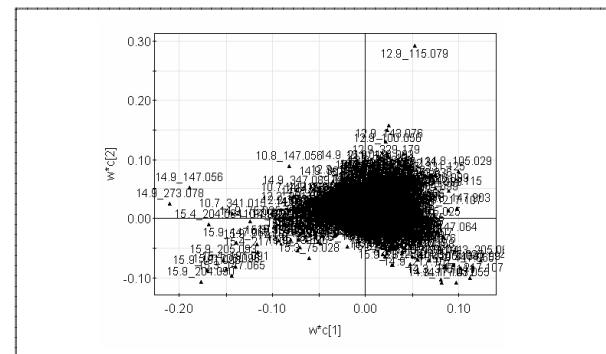


Figure 5.

Application Note

The EI GC/MS data was subjected to peak deconvolution and integration followed by PCA using the MarkerLynx™ Application Manager for MassLynx™ Software, Figure 3. A total of 1282 markers were detected in the EI GC/MS data set. The subsequent PCA scores plot of component 1 versus component 2 is displayed in Figure 4. We can see from this data that the four animal groups (control week 6, Zucker week 6, control week 20, and Zucker week 20) are easily separated into distinct regions. This result demonstrates that GC/MS can be used in metabolomics applications to determine the differences between biological and temporal groups. The resulting loadings plot from the multivariate analysis of the EI GC/MS data is given in Figure 5. In this example, each analyte is labeled with its retention time and m/z value. The analytes furthest away from the origin of the plot are those which contribute most strongly to the variance observed in the data set. One such ion is m/z = 457. This ion was subjected to elemental composition analysis using the exact mass value of 457.1943, giving a result of

$C_{17}H_{37}N_4O_3Si_4$ with a mass error of 3.3 ppm. This data was then submitted to a NIST library search which resulted in a hit for TMS-derivatized uric acid (match score of 792, probability of 82%) shown in Figure 6. The agreement between the exact mass measurement and library search results is excellent and is supported by the CI data.

CONCLUSION

We have demonstrated the utility of GC/MS(TOF) for metabolomic studies of biological fluids following sample derivatization. After data deconvolution and integration, the resulting statistical data indicates differences as a result of disease or toxic insult. One of the significant advantages of GC/MS is that the ions of interest identified in the loadings plot can be subsequently searched using the NIST database for analyte identification.

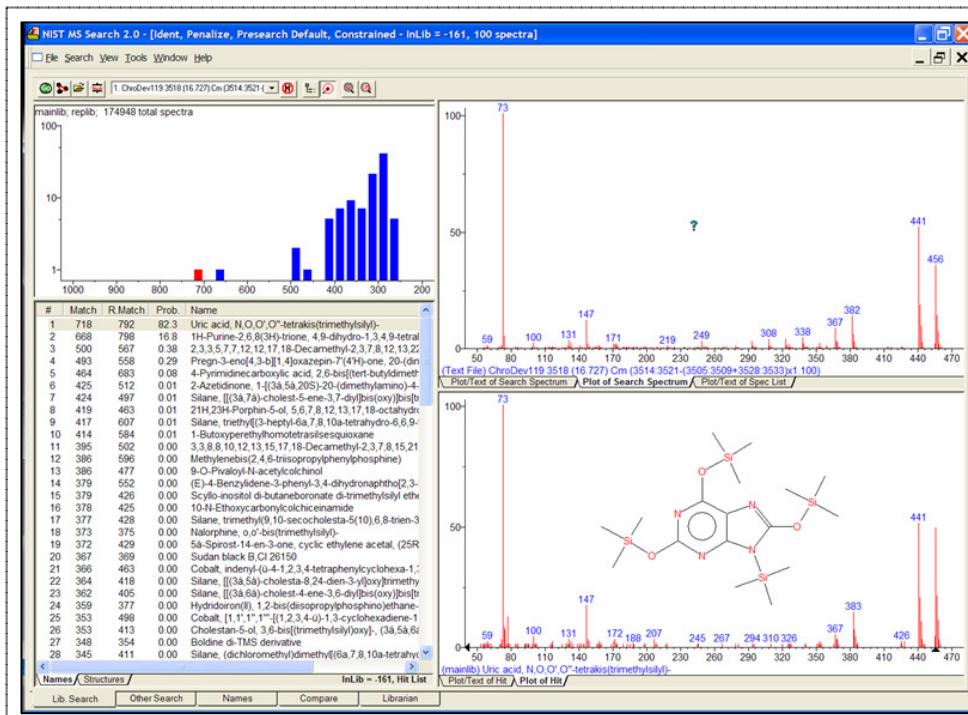


Figure 6.

AUSTRIA AND EXPORT (CENTRAL EUROPE, CIS,
MIDDLE EAST, INDIA AND INDIA SUBCONTINENT)
43 1 877 18 07

AUSTRALIA 61 2 9933 1777

BELGIUM AND LUXEMBOURG
32 2 726 1000

BRAZIL 55 11 5543 7788

CANADA 800 252 4752 X2205

CZECH REPUBLIC 420 2 617 11384

DENMARK 45 46 59 8080

FINLAND 358 09 506 4140

FRANCE 33 1 3048 7200

GERMANY 49 6196 400600

HONG KONG 852 29 64 1800

HUNGARY 36 1 350 5086

INDIA 91 80 837 1900

IRELAND 353 1 448 1500

ITALY 39 02 27 4211

JAPAN 81 3 3471 7191

KOREA 82 2 3284 1300

MEXICO 52 55 5524 7636

THE NETHERLANDS 31 76 508 7200

NORWAY 47 6 384 6050

PEOPLE'S REPUBLIC OF CHINA
86 10 8451 8918

POLAND 48 22 833 4400

PUERTO RICO 787 747 8445

RUSSIA/CIS 7 095 931 9193

SINGAPORE 65 6278 7997

SPAIN 34 93 600 9300

SWEDEN 46 8 555 11 500

SWITZERLAND 41 62 889 2030

TAIWAN 886 2 2543 1898

UK 44 208 238 6100

U.S.A. AND ALL OTHER COUNTRIES:
WATERS CORPORATION
34 Maple St.
Milford, MA 01757 U.S.A.
T: 508 478 2000
F: 508 872 1990
www.waters.com

Waters

WATERS CORPORATION
34 Maple St.
Milford, MA 01757 U.S.A.
T: 508 478 2000
F: 508 872 1990
www.waters.com

For Complete Confidence

Waters, Micromass, GCT, MassLynx, and MarkerLynx are trademarks of Waters Corporation.
All other trademarks are property of their respective owners.
©2004 Waters Corporation. Produced in the U.S.A.
Dec. 2004 720001075EN KJ-PDF

