

ACCURACY AND PRECISION OF A UPLC™/MS/MS SYSTEM FOR THE ANALYSIS OF CAFFEINE IN HUMAN PLASMA

Ed Sprake and Iain Gibb

Waters Corporation, Manchester, UK

INTRODUCTION

Caffeine is a very common analyte within bioanalytical research, due its use as a CYP450 marker and its common application in co-dosing studies. In this technical note, we show the analysis of caffeine in protein precipitated human plasma, performed using the Waters® ACQUITY Ultra Performance LC™ System coupled with the Waters Micromass® Quattro Premier™ Mass Spectrometer. In this work, we have used an external calibration method to highlight the excellent accuracy and precision of this system configuration.

EXPERIMENTAL

Caffeine spiking solutions were made up at 14 concentrations between 45 and 7500 ng/mL in water as shown in Table 1.

First, 100 µL of spiking solution was added to 200 µL of human plasma. Next, 1.0 mL of acetonitrile was added to perform the protein precipitation, and the resulting mixture was spun down at 13,000 rpm for 5 minutes. Finally, 200 µL of the supernatant was diluted into 1.0 mL of the initial mobile phase.

Calibration standards were injected in duplicate, 6 separate quality control (QC) samples were prepared at each of the QC concentration points. The QC samples were injected once each.

LC Conditions

LC System: ACQUITY UPLC™ system
Column: ACQUITY BEH C₁₈
2.1 x 50 mm, 1.7 µm
Eluent: 85:15 water/acetonitrile (isocratic)
Run Time: 0.6 min
Injection Volume: 5 µL

MS Conditions

MS System: Quattro Premier tandem quadrupole mass spectrometer
Ion Mode: Electrospray positive
Capillary Voltage: 3.00 kV
Cone Voltage: 25 V
Detection Mode: MRM (195.1>137.9 Da)
Collision Energy: 20 V
Dwell Time: 0.1 sec
Collision Gas: Argon (3.4x10⁻³ mbar)

Spike Conc. (ng/mL)	Actual Conc. in Plasma (ng/mL)	Sample Type
45	30	Standard
105	70	
300	200	
645	430	
1125	750	
1800	1200	
2700	1800	
3600	2400	
4650	3100	
6000	4000	
7500	5000	QC
45	30	
105	70	
150	100	
450	300	
3000	2000	
7500	5000	

Table 1. Caffeine spiking solutions.

RESULTS AND DISCUSSION

A full validation batch was run on each of 3 separate days. A typical calibration curve is shown in Figure 1. The calibration curve was plotted using a linear fit with a $1/x^2$ weighting, and gave a coefficient of correlation of >0.998.

Shown in Tables 2 and 3 are the inter-batch statistics for the standard and QC samples. The statistics for the standard injections are based on 2 replicate injections of the 11 calibration points for each of the 3 inter-day batches. All calibration points show <7% CV with accuracy values between 97.9% – 102.8% (Table 2). The statistics for the QC injections are based on single injections of 6 individually spiked QC solutions at each concentration for each of the 3 inter-day batches. All QC levels show <6.4% CV with accuracy values between 98.1% – 100.9% (Table 3).

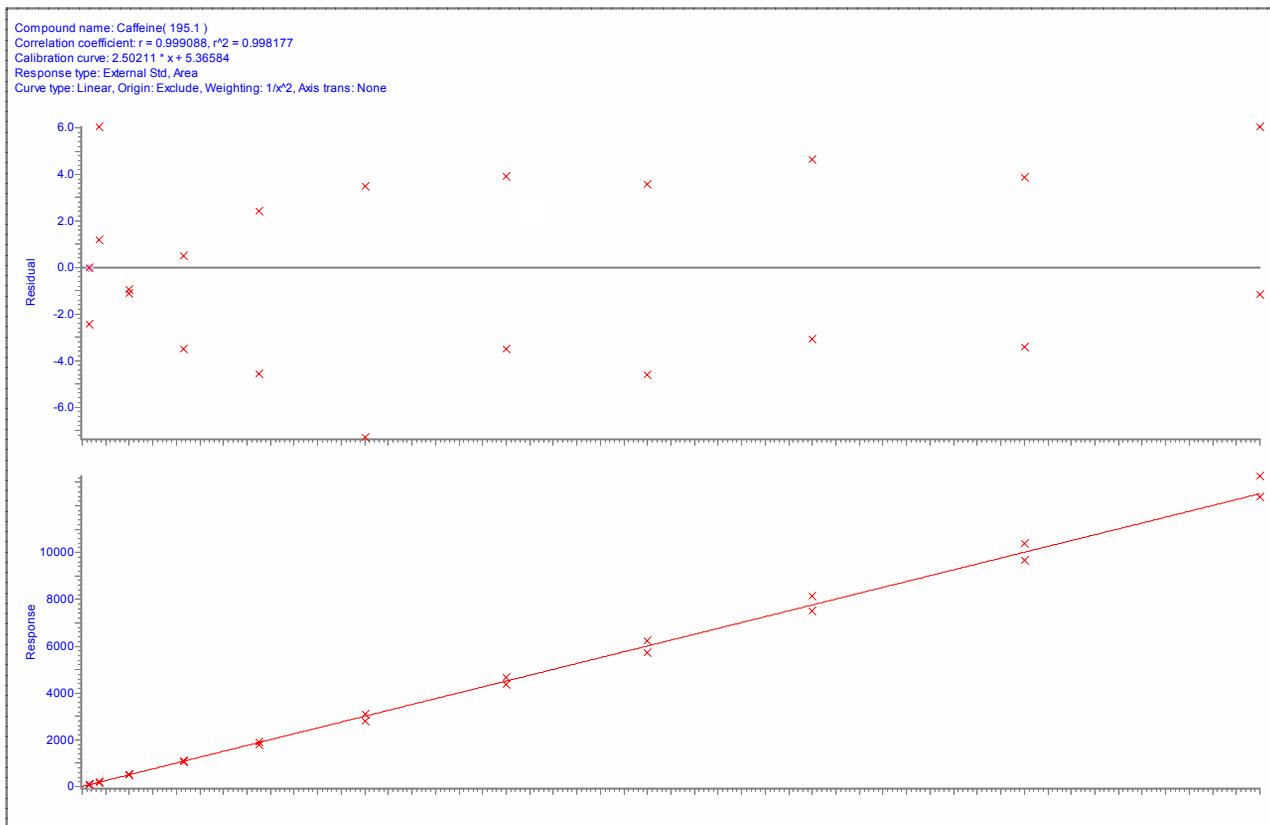


Figure 1. Calibration curve of caffeine in human plasma showing a value for R^2 of >0.998.

Conc. of Caffeine (ng/mL)	Mean (n=6)	SD (n=6)	%CV (n=6)	%Accuracy (n=6)
30	30.0	2.1	6.96	99.9
70	71.1	1.8	2.50	101.5
200	195.7	12.6	6.46	97.9
430	421.6	23.8	5.64	98.0
750	740.5	33.0	4.46	98.7
1200	1187.4	58.5	4.92	98.9
1800	1798.2	83.1	4.62	99.9
2400	2393.8	101.8	4.25	99.7
3100	3133.4	110.9	3.54	101.1
4000	4060.5	162.7	4.01	101.5
5000	5142.1	184.9	3.60	102.8

Table 2. Inter-batch statistics for the calibration standard injections across three inter-day validation batches.

CONCLUSIONS

The ACQUITY UPLC system with the Quattro Premier tandem quadrupole mass spectrometer have been combined to yield an unmatched platform for quantitative LC/MS/MS. The results demonstrate the excellent accuracy and precision that can be achieved for compounds in complex biological matrices, such as human plasma. In this case, the required criteria were easily met without the use of an internal standard. With the enhanced chromatographic capabilities of UPLC, this analysis was achieved with a run time of 0.6 minutes. This shows that accuracy and precision can be achieved with very fast run times, giving the potential to significantly increase throughput as compared to an HPLC-based system, yet still maintain the necessary selectivity and resolution for good quality data. The target LLOQ of 30 ng/mL is much higher than could potentially be achieved with UPLC/MS/MS, but was determined to be sufficient for the dosage levels to be detected.

WATERS CORPORATION
34 Maple St.
Milford, MA 01757 U.S.A.
T: 508 478 2000
F: 508 872 1990
www.waters.com

Conc. of Caffeine (ng/mL)	Mean (n=18)	SD (n=18)	%CV (n=18)	%Accuracy (n=18)
30	29.6	1.87	6.32	98.6
70	68.9	3.66	5.31	98.4
100	99.7	3.67	3.69	99.7
300	294.3	11.2	3.80	98.1
2000	2018.0	113.2	5.61	100.9
5000	5031.4	136.3	2.71	100.6

Table 3. Inter-batch statistics for the QC injections across three inter-day validation batches.

Waters

For Complete Confidence

Waters, Micromass, ACQUITY UPLC, UPLC, Ultra Performance LC, and Quattro Premier are trademarks of Waters Corporation.

All other trademarks are the property of their respective owners.

©2005 Waters Corporation Produced in the U.S.A. Nov. 2005 720001413EN KJ-PDF

