

ENHANCED DETERMINATION OF THE METABOLIC FATE OF DRUGS USING UPLC™/oa-TOF MS

Mike McCullagh, Jose Castro-Perez, and Lisa Calton
Waters Corporation, Manchester, UK

INTRODUCTION

The application area of drug metabolism analysis is complex and extremely challenging. LC/MS is well-suited for this, due to its inherent sensitivity and selectivity. It has revolutionized many areas of the drug discovery process, in particular lead optimization, where tandem mass spectrometry is widely used to generate ADME information for new compounds. Quantitative data is usually provided by MRM experiments using tandem quadrupole MS to determine rate of metabolism, but does not provide qualitative information on the sited mechanism of metabolism. The prevalent requirement to identify metabolites has led to a significant increase in the use of exact mass measurements. LC coupled to orthogonal acceleration-time of flight (oa-TOF) mass spectrometers addresses this need by providing sensitive and accurate qualitative and quantitative information from one analysis.

The dynamic range of LC-TOF instruments has been improved, removing errors in quantification and exact mass measurement at high metabolite concentrations.

Combining the Waters® Micromass® Q-ToF Premier™ Mass Spectrometer with the Waters ACQUITY Ultra Performance LC™ System, which provides faster analysis and improved chromatographic resolution, we can generate higher quality data to permit automated data processing and a unique mass defect to confidently distinguish drug-related metabolites from matrix components.

In this work, we present data using a UPLC™/oa-TOF MS platform with extended dynamic range capability that broadens its utility for metabolism studies. Rat microsomal incubations of 5 drug candidates were analyzed: Verapamil, Clozapine, Glyburide, Dextromethorphan and Ketoconazole (due to the extensive data produced, only verapamil and dextromethorphan are shown). The metabolites were automatically identified using the Metabolynx™ Application Manager for MassLynx™ Software, a post-acquisition data processing package in which expected and unexpected metabolites can be identified by comparing control and analyte samples. Metabolynx has been enhanced further to use the specificity of exact mass measurement and apply mass defect filters. The

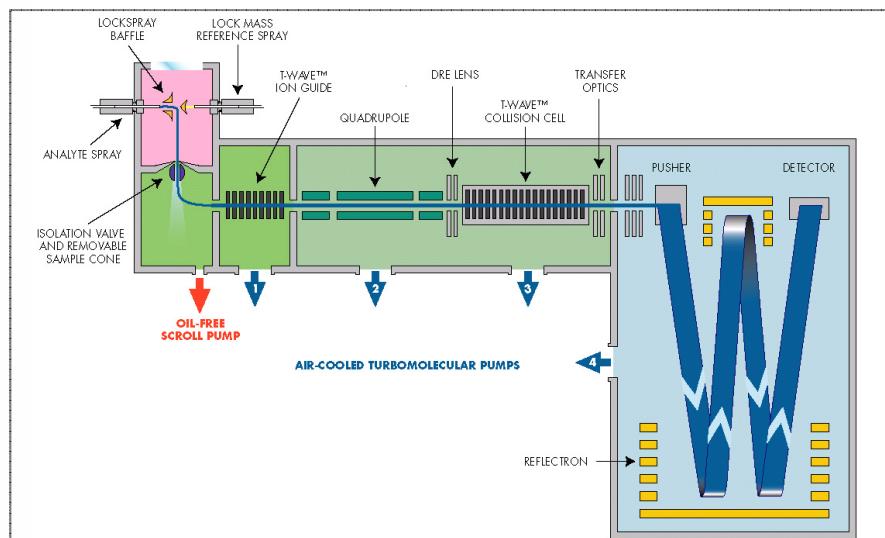


Figure 1. Q-ToF Premier mass spectrometer schematic.

mass defect filter utilizes the exact mass and mass deficiency of the parent drug compound to rapidly identify only drug-related metabolites.

EXPERIMENTAL

MS Conditions

MS System:	Q-ToF Premier oa-TOF mass spectrometer
Ionization Mode:	ESI ⁺ at 3 kV
Cone Voltage:	40 V
Reference Mass:	Leucine enkephalin, [M+H] ⁺ = 556.2771
Acquisition Details:	100 – 1000 m/z, 0.2 sec/spectrum 0.1 sec inter-acquisition delay
Resolution:	>8000 FWHM (V mode)

UPLC Conditions

LC System:	ACQUITY UPLC system
Column:	ACQUITY UPLC BEH C ₁₈ 2.1 X 100 mm, 1.7 µm
Column Temp:	45 °C
Flow:	0.6 mL/min
Mobile Phase:	A: H ₂ O (0.1% HCOOH) B: MeCN
Gradient:	0.0 min 2% B, 8.0 min 98% B, 10.0 min 98% B, 10.1 min 2% B, 13.0 min 2% B

HPLC Conditions

LC System:	Waters Alliance® HT system
Column:	Symmetry® C ₁₈ 2.1 X 100 mm, 3.5 µm
Column Temp:	35 °C
Flow:	0.3 mL/min
Mobile Phase:	A: H ₂ O (0.1 % Formic Acid) B: MeCN
Gradient:	0.0 – 0.5 min 2% B, 0.5 – 20.0 min 80% B, 20.0 – 30.0 min 80% B, 30.0 – 30.5 min 2% B, 30.5 – 35.0 min 2% B

Microsomal Incubation Conditions

Microsomes:	Human microsomes/S9
Water Bath Temp:	37 °C
Buffers:	50 – 100 mM Tris or K Phosphate @ pH 7.4
NADPH Regenerating System (NRS) in 2% NaHCO ₃ (Sigma S-5761) Solution:	0.5 mg/mL b-NADP (Sigma N-0505), 2.0 mg/mL Glucose-6-phosphate (Sigma G-7879), 1.5 units/mL Glucose-6-phosphate dehydrogenase (Sigma G-7877)
Quench:	Equal volume MeCN
Drug Concentration:	100 µL 0.1 mg/mL, final quench volume 2 mL

RESULTS

In Figure 2, the UPLC/TOF extracted mass chromatograms for metabolites formed for the *in vitro* incubation of dextromethorphan are shown at T = 15 min, T = 30 min, and T = 60 min time points. Presented in Figure 3 are the exact mass spectra for the demethylated metabolite formed for the *in vitro* incubation of dextromethorphan at each of the time points described above. Illustrated in Figure 4 are the exact mass spectra for the parent drug at each of the time points. In Figure 5, the exact mass spectra are shown for the hydroxylated metabolite formed for the *in vitro* incubation of dextromethorphan at each of the time points. The HPLC/TOF/MS extracted mass chromatograms for metabolites formed for the *in vitro* incubation of verapamil are presented in Figure 6. The corresponding UPLC/TOF/MS extracted mass chromatograms are summarized in Figure 7, with the un-metabolized parent (labelled "A"). The metabolites shown are identified in Table 1 (labeled "B" through "J").

Figure	Metabolite Identified
7b.	cleaved N-de-alkylation/hydroxylation
7c.	cleaved/de-methylation
7d.	cleaved
7e.	glucuronidation/de-ethylation
7f.	de-ethylated
7g.	glucuronidation/de-methylation
7h.	hydroxylation
7i.	de-alkylation/hydroxylation
7j.	de-methylation

Table 1. A list of the metabolites identified in Figure 7.

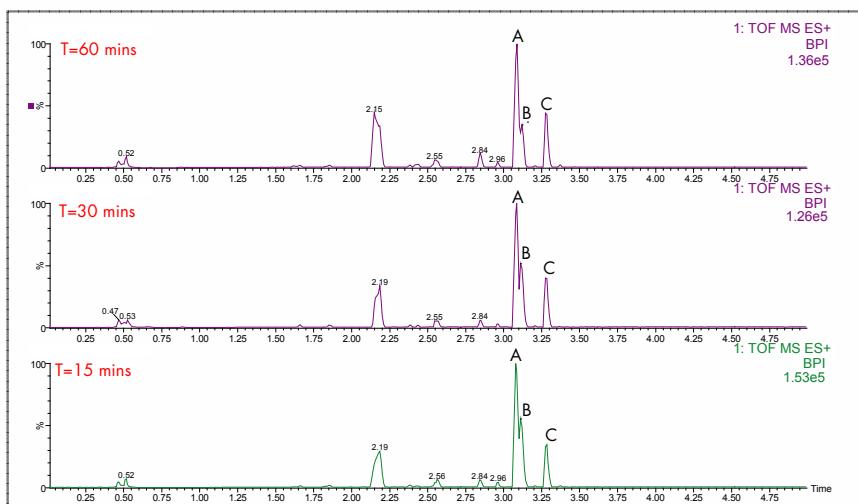


Figure 2. UPLC/TOF extracted mass chromatograms for metabolites formed for the *in vitro* incubation of dextromethorphan at time points 15 min, 30 min, and 60 min.

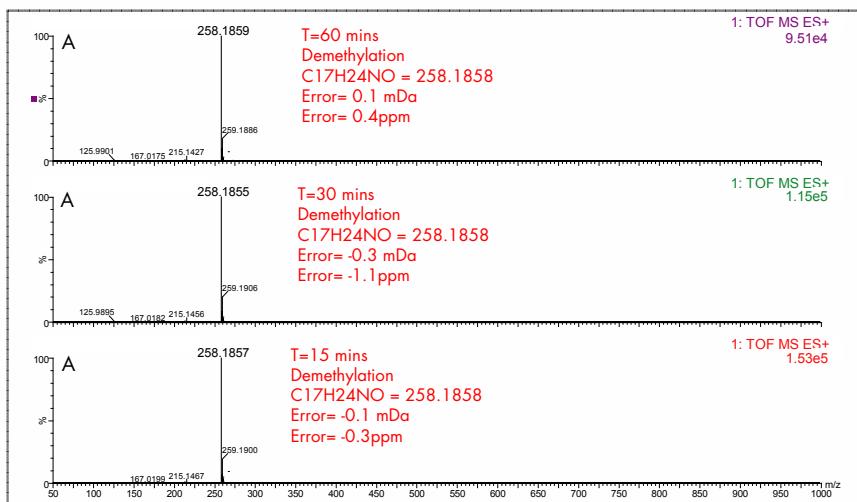


Figure 3. Exact mass spectra for the demethylated metabolite formed for the *in vitro* incubation of dextromethorphan at time points 15 min, 30 min, and 60 min.

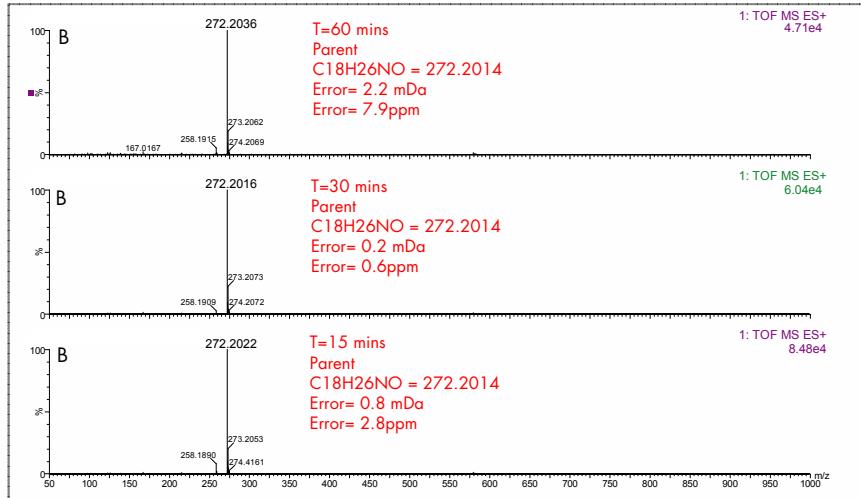


Figure 4. Exact mass spectra for the *in vitro* incubation of the parent dextromethorphan drug at time points 15 min, 30 min, and 60 min.

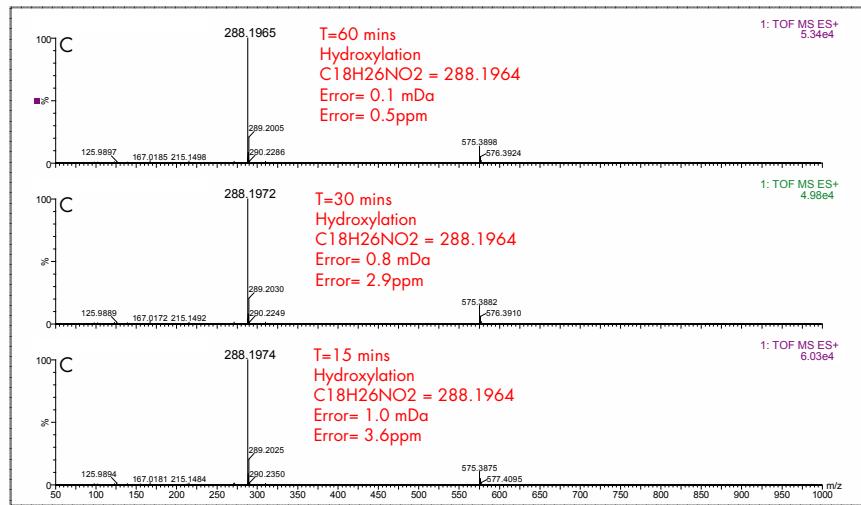


Figure 5. Exact mass spectra for the hydroxylated metabolite formed for the *in vitro* incubation of dextromethorphan at time points 15 min, 30 min, and 60 min.

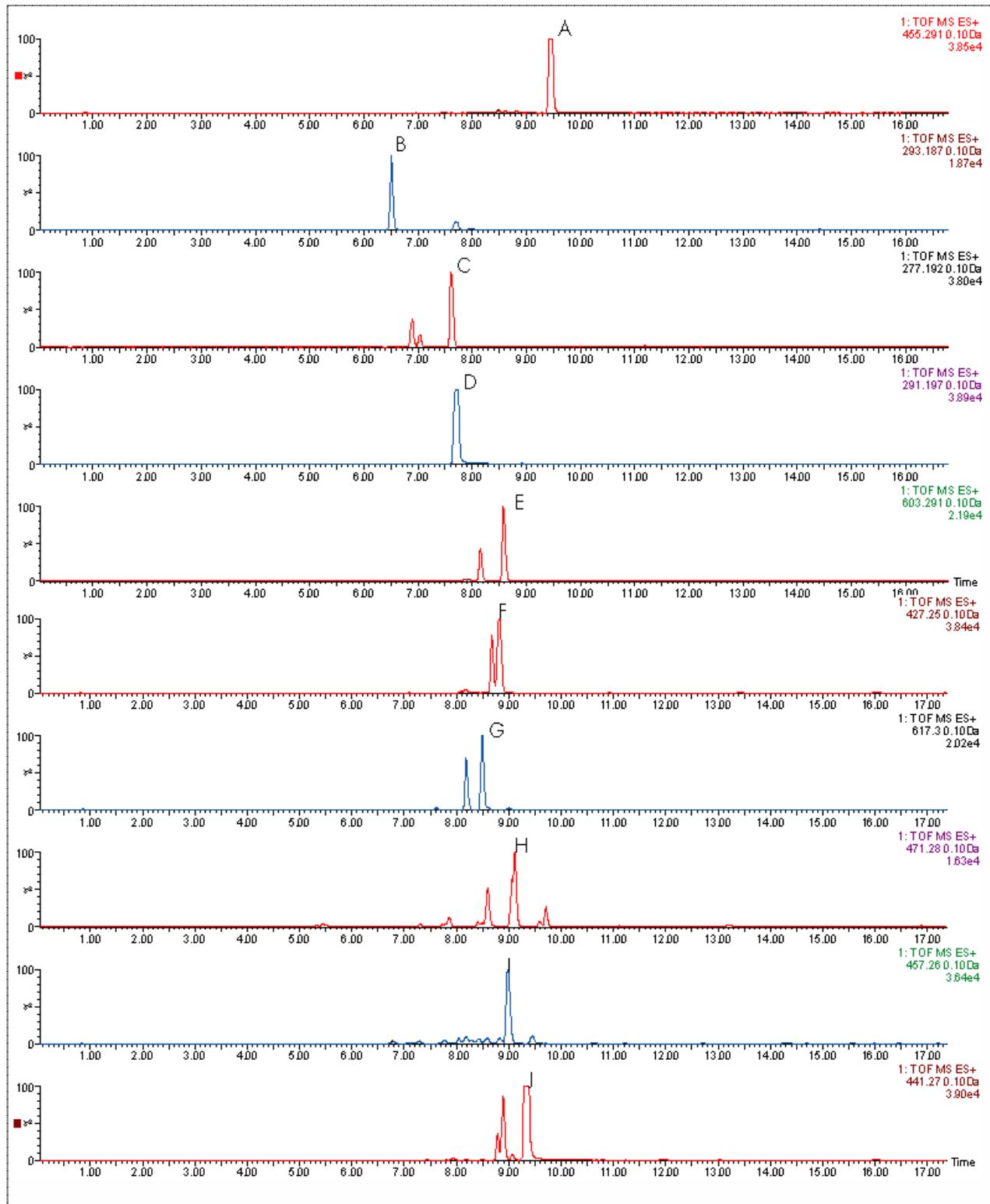


Figure 6. HPLC/oa-TOF MS extracted mass chromatograms for metabolites formed for the *in vitro* incubation of verapamil.

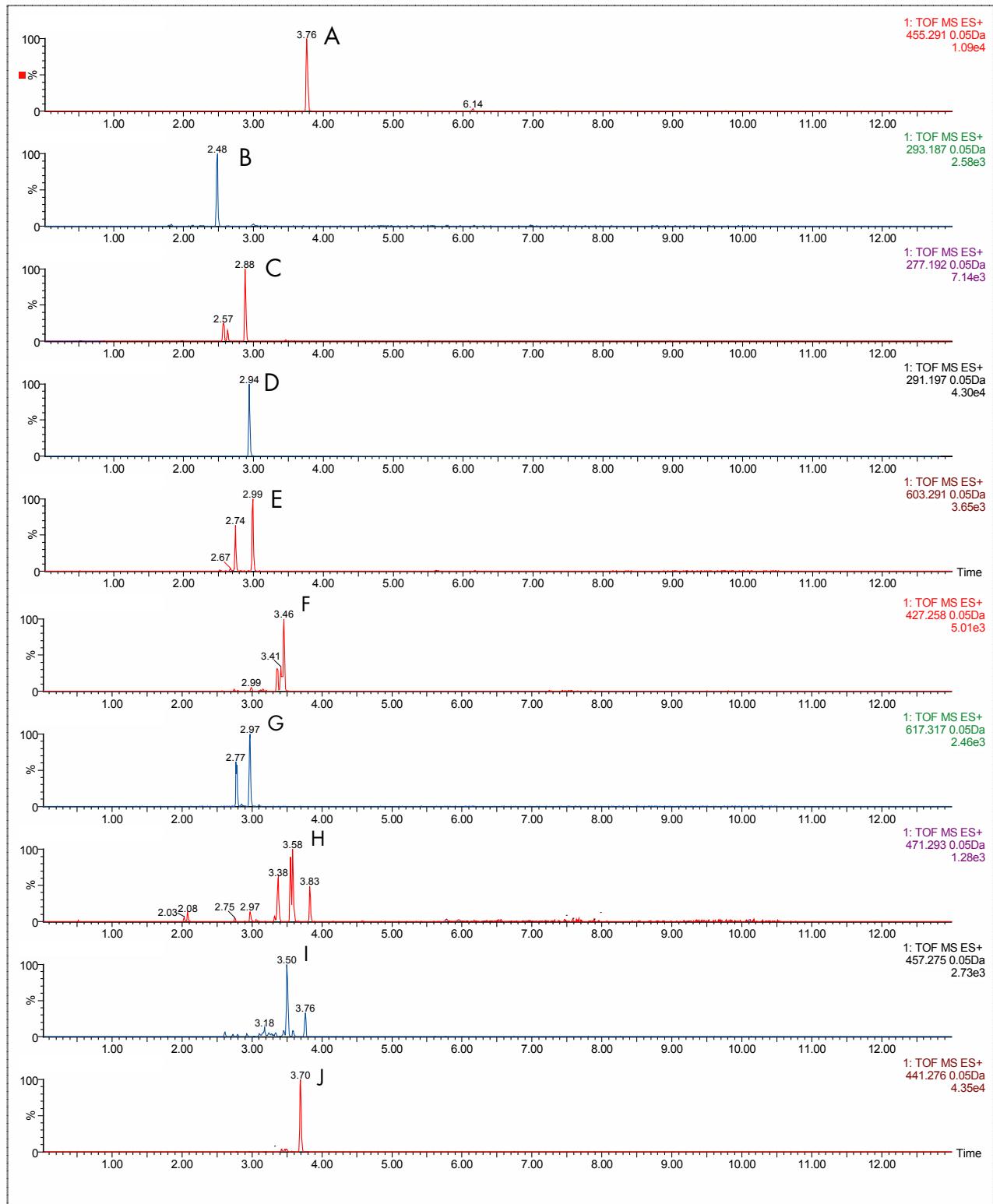


Figure 7. UPLC/oa-TOF MS extracted mass chromatograms for metabolites formed for the *in vitro* incubation of verapamil.

The reduced data set produced by MetaboLynx is displayed in a browser format in Figure 8, where the results of application of the post-processing mass defect data filter are illustrated.

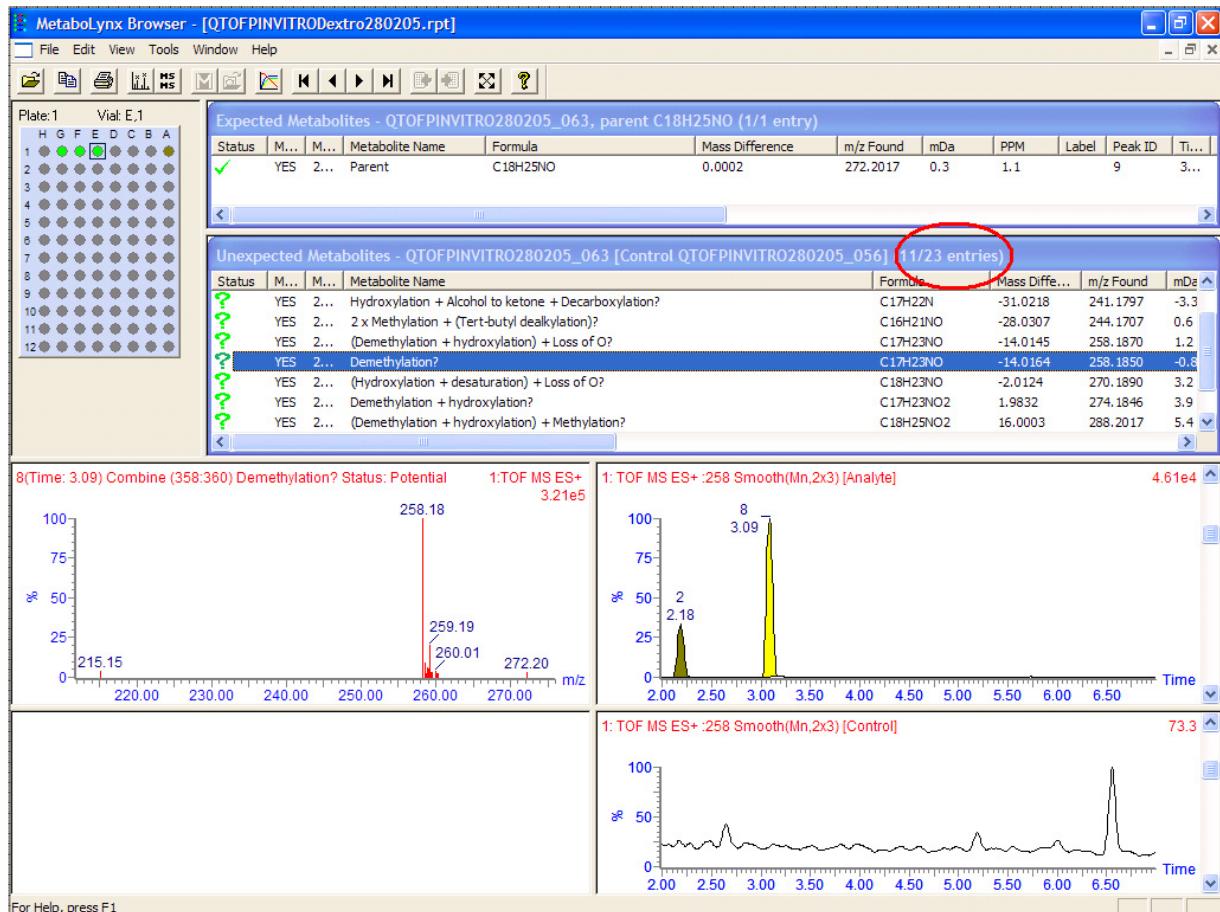


Figure 8. The MetaboLynx report browser illustrating application of post-processing mass defect data filter.

DISCUSSION

The parent drugs verapamil and dextromethorphan were incubated *in vitro* using human S9 microsomes. These incubations were utilized to test both hardware and software enhancements for determining the metabolic fate of drugs. The schematic of a Q-Tof Premier in Figure 1 shows the pDRE™ lens, which enhances the analytical applicability of oa-TOF MS. This feature has enabled extended dynamic range up to 4 orders of magnitude, providing both qualitative and quantitative capabilities with good exact mass measurements for highly concentrated peaks. Previously, at high ion counts, a shift to lower mass would occur, giving incorrect mass measurement and less confidence in elemental composition determination. Selection of spectra at low ion counts would provide excellent mass measurement and high confidence in the resultant data, however, this could be a laborious task. Even at high ion counts, pDRE allows either the peak top, or the whole peak to be averaged to generate an exact mass spectrum. This improves the ease with which data can be generated and processed. Also, since no detector saturation is taking place, a true profile of the route and rate of metabolism will be obtained. The impact of dynamic range enhancement can be seen in Figure 2 where the UPLC/TOF BPI chromatograms for metabolites formed, for the *in vitro* incubation of

dextromethorphan are shown at time points 15 min, 30 min, and 60 min. A true chromatographic profile is observed. The exact mass spectra obtained for the demethylated (labeled "A") and hydroxylated metabolites (labeled "C"), can be seen for 3 time point intensities in Figures 3 and 5, respectively. In each case the spectra comprised of a single spectrum, generated from the top of the peak. Intensities of up to 150,000 counts per ion accumulation are illustrated, where no detector saturation has been observed and mass measurement errors of typically less than 3 ppm have been obtained. The corresponding spectra produced for the remaining parent dextromethorphan are shown in Figure 4. pDRE has enabled exact mass measurement to be obtained routinely over a suitable concentration range for true measurement of metabolite formation and profiling.

As well as improving mass measurement performance, chromatography technology has also been enhanced by using Ultra Performance LC. In Figures 6 and 7, the HPLC and UPLC/TOF extracted mass chromatograms for metabolites formed for the *in vitro* incubation of verapamil are presented. The HPLC data was acquired as part of a previous study.¹ It is shown that not only is analysis time reduced by greater than 50%, but chromatographic resolution is also improved. Metabolite isomers can now be chromatographically resolved, reducing the risk of "missed" metabolites. Improved chromatographic resolution also adds further opportunity to obtain higher quality MS data, reducing the need for multiple simultaneous MS experiments.

An illustration of the Metabolynx report browser is contained in Figure 8. The processing of metabolism data is usually a significant bottleneck in metabolite identification. Using Metabolynx to automatically process the data accelerates this step. The reduced data are displayed on a results browser, which allows rapid interpretation of the data. In addition the data can be greatly simplified by application of the post-processing mass defect data filter. The mass defect

property of the parent drug will be maintained within a narrow mass range window, for both phase I and phase II biotransformations. The ability to acquire enhanced quality accurate mass spectra provides the opportunity to exploit mass sufficiency and deficiency. Application of a +/- 7 mDa mass defect window has enabled the number of unexpected metabolites found to be reduced to 11 from 23. This post-processing step enabled immediate removal of 12 non drug matrix-related unexpected metabolites to be removed.

CONCLUSION

- Analysis time reduced by at least 50% using a generic method UPLC method.
- Improved chromatographic resolution has been obtained routinely.
- pDRE extends the dynamic range for which accurate mass measurement can be made to 4 orders, enabling specific routine route and rate determination.
- Mass measurement errors of < 3 ppm have been obtained routinely at high ion intensities.
- The highly specific nature of exact mass measurement enables elemental compositions to rapidly be determined and therefore, absolute confidence in producing the correct identification of metabolites formed.
- Application of a post-processing mass defect data filter enables rapid exclusion of non drug-related unexpected metabolites.

REFERENCES

1. Mike McCullagh, Jose Castro-Perez, Andrew Baker, and Jennifer Granger. Simultaneous determination of route and rate metabolism using LC-TOF. Poster presented at the ISSX Drug Metabolism Meeting in Vancouver, B.C., August 29 – September 1, 2004. Waters Literature Code 720000972EN.

Sales Offices:

Austria 43 1 877 18 07

Australia 61 2 9933 1777

Belgium and Luxembourg

32 2 726 1000

Brazil 55 11 5543 7788

Canada 800 252 4752

China 86 10 8451 8918

Czech Republic 420 2 617 11384

Denmark 45 46 59 8080

Finland 358 9 506 4140

France 33 1 3048 7200

Germany 49 6196 400600

Hong Kong 852 29 64 1800

Hungary 36 1 350 5086

India 91 80 2837 1900

Ireland 353 1 448 1500

Italy 39 02 27 4211

Japan 81 3 3471 7191

Korea 82 2 820 2700

Mexico 52 55 5200 1860

The Netherlands 31 76 508 7200

Norway 47 6 384 6050

Poland 48 22 833 4400

Puerto Rico 787 747 8445

Russia/CIS 7 095 931 9193

Singapore 65 6278 7997

Spain 34 93 600 9300

Sweden 46 8 555 11 500

Switzerland 41 62 889 2030

Taiwan 886 2 2543 1898

UK 44 208 238 6100

US 1 800 252 4752

WATERS CORPORATION

34 Maple St.

Milford, MA 01757 U.S.A.

T: 508 478 2000

F: 508 872 1990

www.waters.com

Waters

For Complete Confidence

Waters, MassLynx, Micromass, Metabolynx, Ultra Performance LC, ACQUITY UPLC, UPLC,

Q-ToF Premier, Symmetry, pDRE, and Alliance are trademarks of Waters Corporation.

All other trademarks are the property of their respective owners.

©2005 Waters Corporation Produced in the U.S.A. Dec. 2005 720001419EN KJ-PDF

