

Andrew J. Aubin and Katherine Hynes
Waters Corporation, Milford, MA, USA

INTRODUCTION

Riboflavin (vitamin B2) is a water soluble B-complex vitamin important in human and animal nutrition. Riboflavin metabolizes into two coenzymes, Flavin Adenine Dinucleotide (FAD) and Flavin Mononucleotide (FMN), and also undergoes photodegradation primarily to 7,8-Dimethylalloxazine (Lumichrome). Previously, HPLC separations of these related compounds have been done in approximately twenty five minutes on 150 mm columns with minimal

resolution and poor peak shapes¹. Separations in less than eight minutes have been achieved with poor resolution between FAD and FMN and indications of co-eluting compounds based on the peak shape². This application note will show the Ultra Performance LC™ (UPLC™) separation of these compounds utilizing photodiode array detection and single quadrupole mass spectrometry (UPLC/PDA/MS) for compound confirmation.

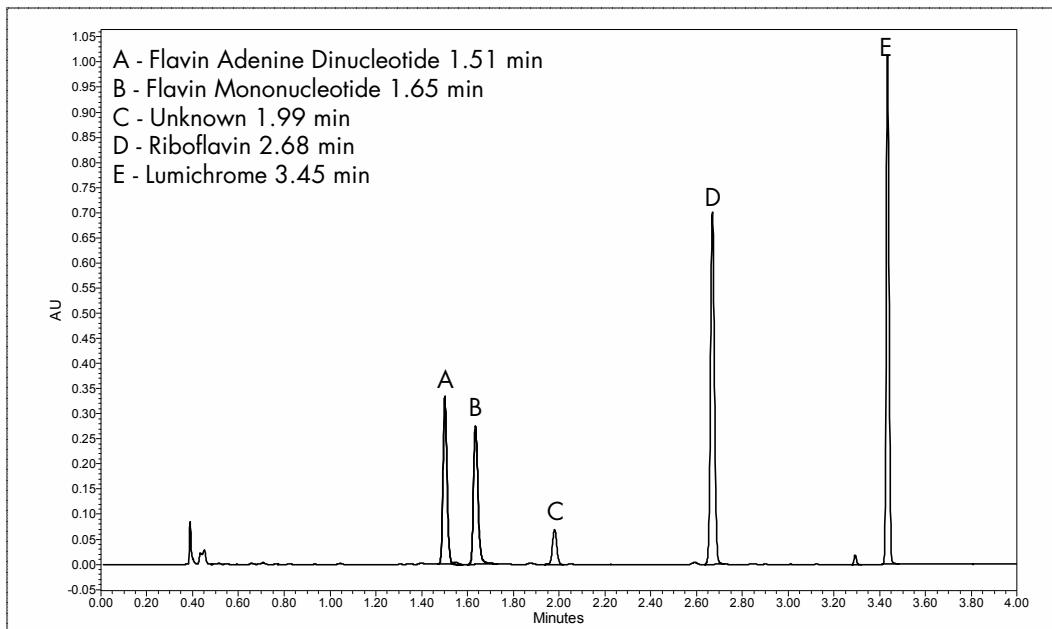


Figure 1. UPLC Separation of Riboflavin, FAD, FMN, and Lumichrome (266 nm).

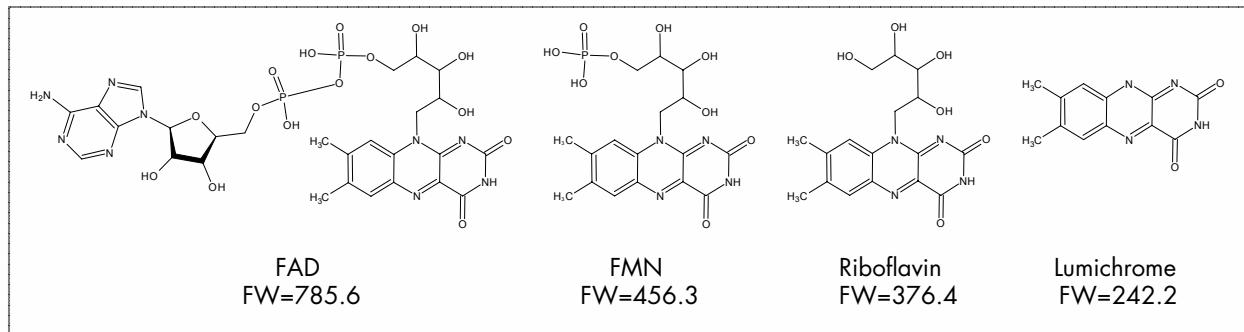


Figure 2. Structures of riboflavin and related compounds.

EXPERIMENTAL

UPLC/PDA/MS Conditions

The Waters® ACQUITY UPLC™ System consisted of a Binary Solvent Manager (BSM) and a Sample Manager (SM) equipped with a 5 μ L loop. Dual detection was employed using an ACQUITY UPLC Photodiode Array Detector in series with a Waters ZQ™ 2000 Mass Detector. For system control, data collection, and processing, Waters Empower™ 2 chromatography data software was employed.

Linear gradient separation (Table 1) of the four related compounds (and other unknown impurities) was achieved on a 2.1 x 100 mm ACQUITY BEH C₁₈ column containing 1.7 μ m particles at a flow rate of 650 μ L/min at 40 °C. Injections of 3 μ L were done using the 'Partial Loop Needle Overfill' mode with the ACQUITY UPLC system. To eliminate carryover, 200 μ L of strong wash (1:1 water/acetonitrile) followed by 600 μ L weak wash (95:5 water/acetonitrile) was used. ZQ 2000 mass detector conditions are described in Table 2.

Time (min)	Profile*	
	%A	%B
0.0	80	20
2.0	60	40
3.0	1	99
3.5	1	99
4.0	80	20

*Where Solvent A = 3.85 g/L Ammonium Acetate in Water and Solvent B = 3.85 g/L Ammonium Acetate in Water/Acetonitrile 7:3.

Table 1. UPLC linear gradient separation conditions.

ES capillary (kV): 3.60	Source Temp (°C): 350
Cone (V): 30	Cone Temp (°C): 120
Extractor (V): 3	Cone Gas Flow (L/Hr): 50
RF Lens (V): 0	Desolvation Flow (L/Hr): 550
LM Resolution: 14	Scan Time: 0.15 s
HM Resolution: 15	Inter-scan Delay: 0.05 s
Multiplier: 650	Scan Range: 150–900 m/z
Probe: ES ⁺	

Table 2. ZQ instrument parameters and tune page parameters.

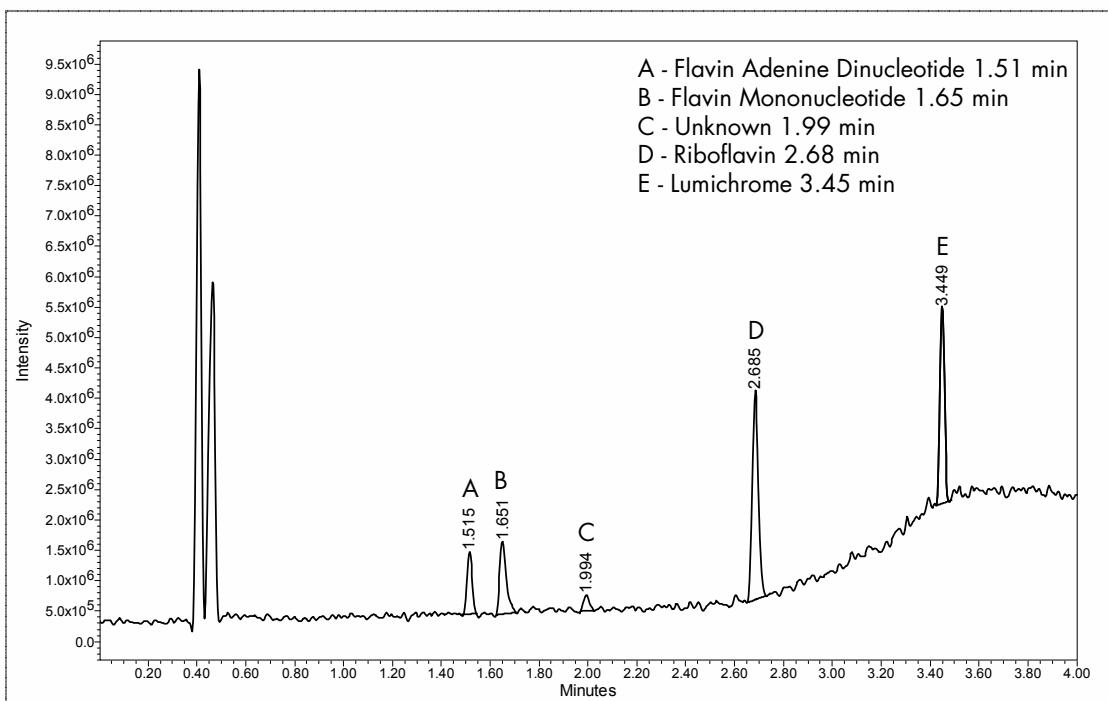


Figure 3. Total ion chromatogram UPLC separation of riboflavin, FAD, FMN and lumichrome.

Sample Preparation

Riboflavin, flavin adenine dinucleotide, flavin mononucleotide, lumichrome, and ammonium acetate were obtained from Sigma-Aldrich Co. Acetonitrile (Optima grade) was obtained from Fisher Scientific. Stock solutions (0.5 mg/mL) of each of the standards were prepared in hot 1M sodium bicarbonate. After cooling, mixed working solutions (0.05 mg/mL of each compound) were prepared in 1M sodium bicarbonate at room temperature. All stock and working solutions were then stored in the dark at 4 °C.

RESULTS AND DISCUSSION

As can be seen in Figures 1 and 3, the four compounds of interest are well separated in under four minutes with UPLC. Examination of the mass spectral data provided positive identification for the four compounds of interest (Figure 4), with the base peak masses matching the molecular weights of the four compounds. In addition to the four identified compounds, another relatively large unknown compound was also resolved from the mixture. This unknown compound (RT=1.98 minutes) had a UV spectrum that was similar to both FMN and riboflavin

as can been seen in Figure 5, indicating that it was likely a related compound. The unknown compound at 1.98 minutes showed a mass of M+H=439.1 (Figure 6).

The generation of ultraviolet (UV) and mass spectra of unknown peaks provides information that can aid the process of identification. In the case of this assay, it was observed that the unknown UV spectrum was very similar to FMN and riboflavin. The unknown mass was consistent with the loss of water from FMN. Its longer retention time is also indicative of this loss. Further identification of this molecule would require the use of more advanced techniques such as time-of-flight mass spectrometry (TOF MS) for exact mass information, and MS/MS for structural confirmation.

Further examination of the PDA data also shows many other well resolved unknown related compounds (Figure 7). No attempt was made to identify these compounds, although the small peak (RT=3.29 min) immediately preceding lumichrome had a UV spectra that was very similar to lumichrome (Figure 8), indicating that it is likely related to lumichrome or is possibly another photodegradation product of riboflavin.

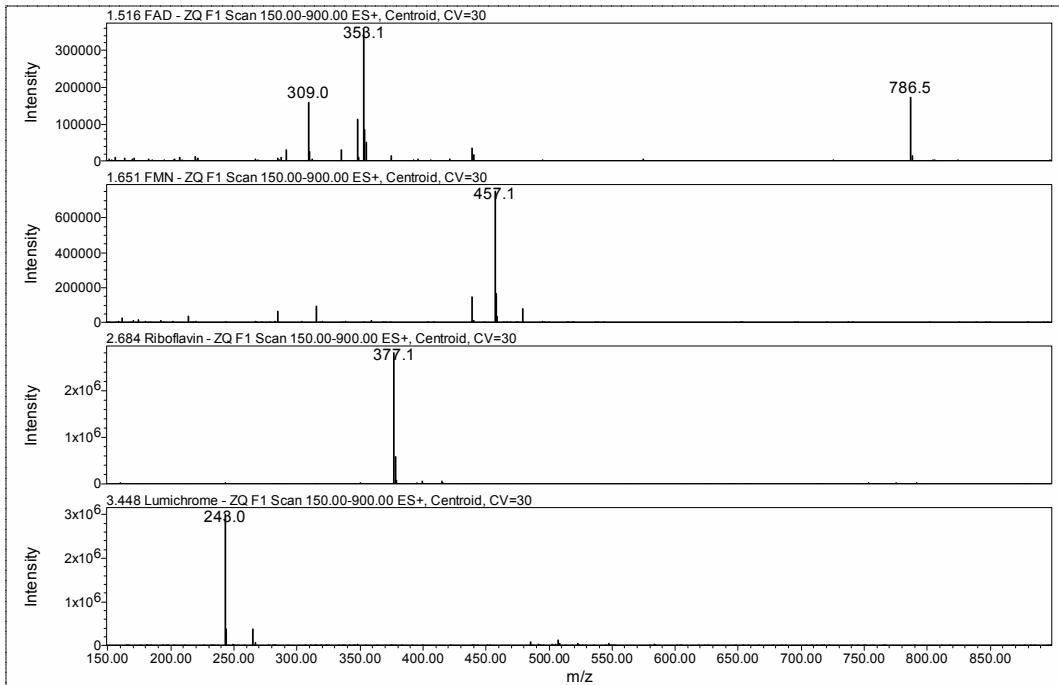


Figure 4. Mass spectra obtained from the ZQ of FAD, FMN, riboflavin, and lumichrome.

Application NOTE

Waters

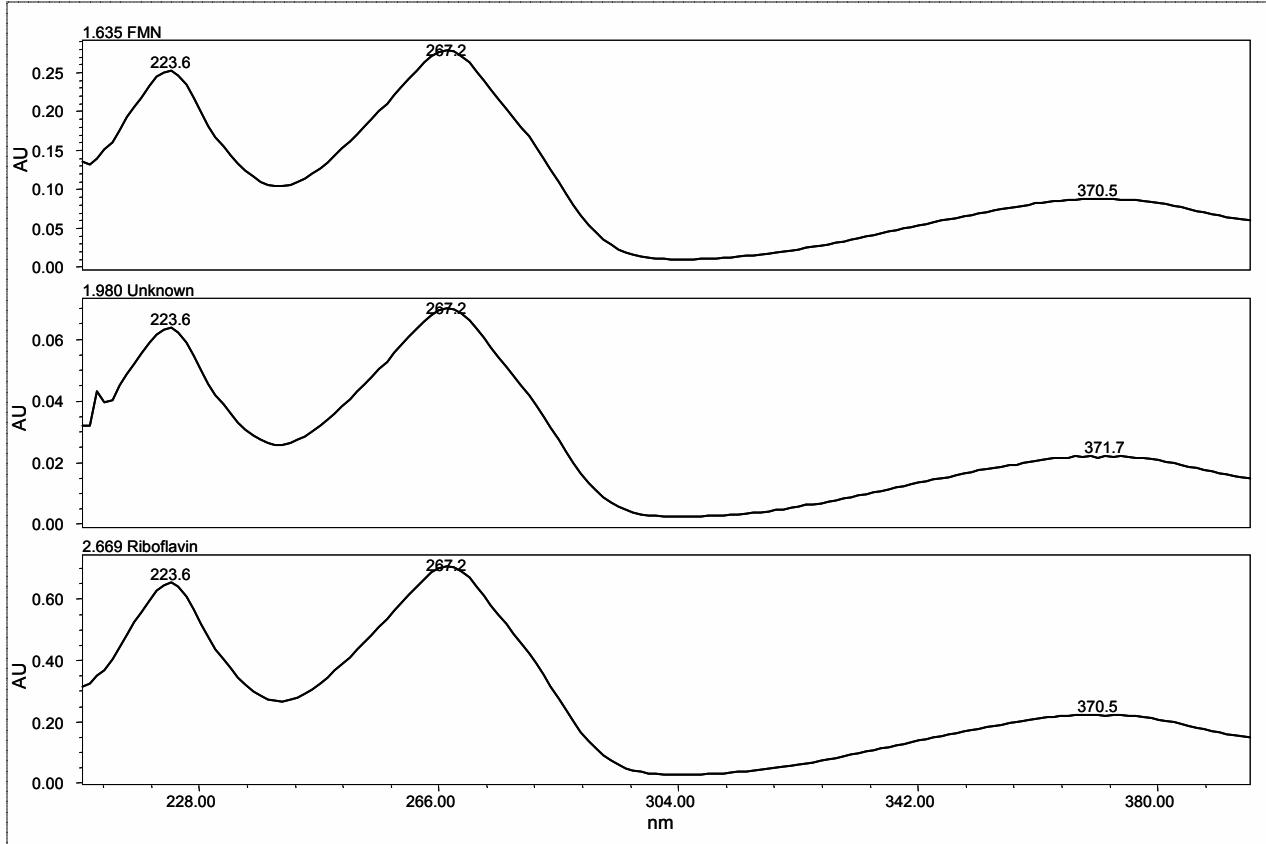


Figure 5. UV spectra of FMN (top), the unknown at 1.98 minutes (middle), and riboflavin (bottom).

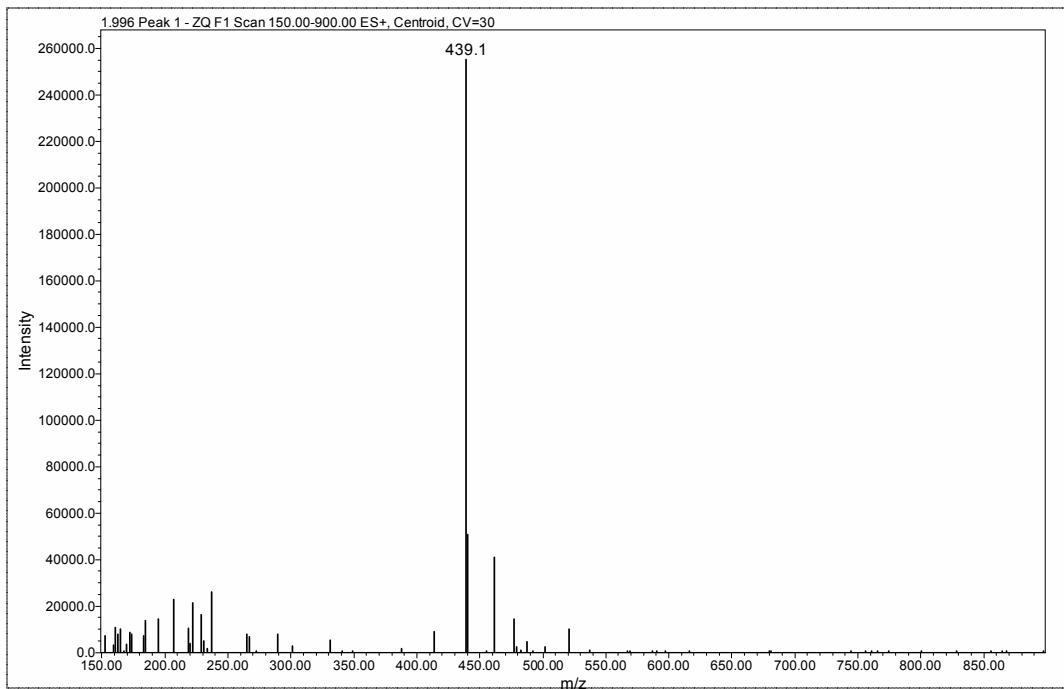


Figure 6. Mass spectrum obtained from the ZQ of the unknown at 1.98 minutes.

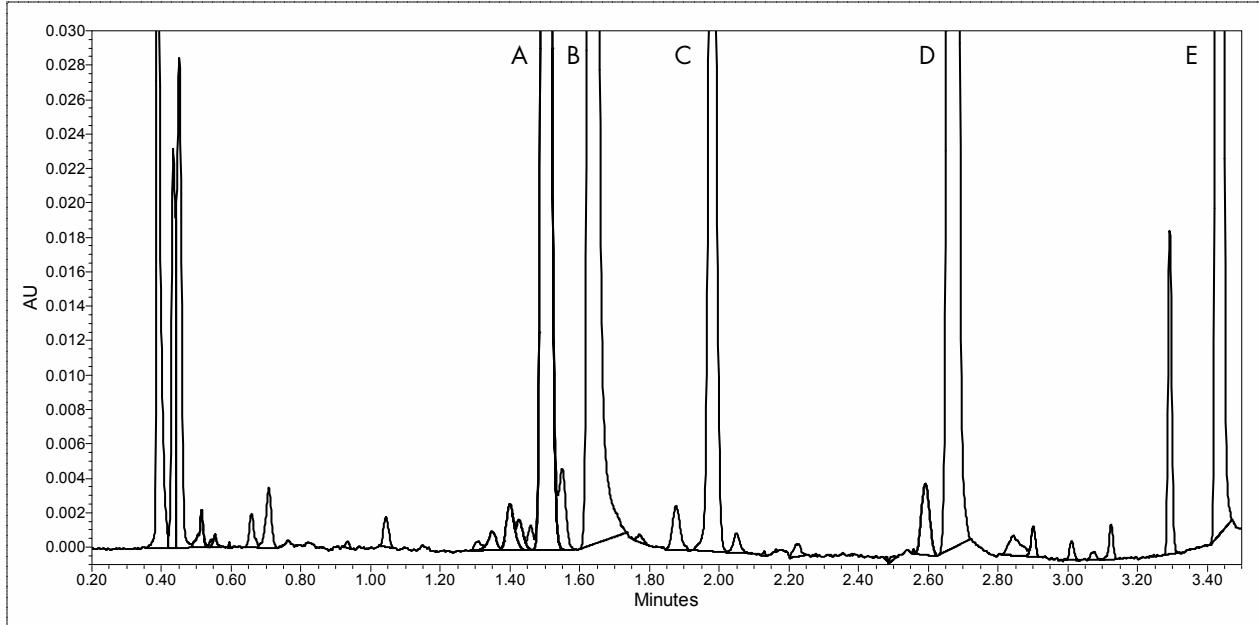


Figure 7. UPLC separation of riboflavin, FAD, FMN and lumichrome (266 nm zoom view).

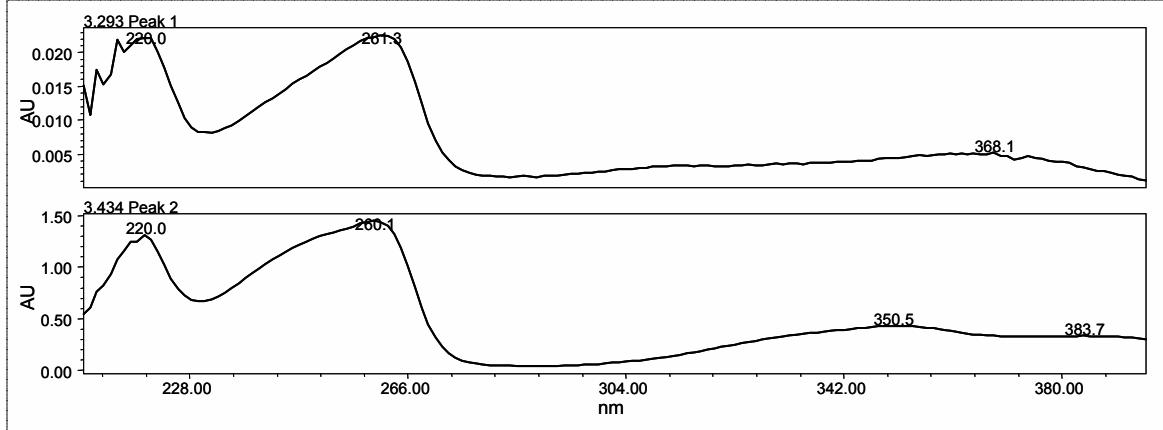


Figure 8. UV spectra of the unknown at 3.29 minutes (top), and lumichrome (bottom).

CONCLUSIONS

The separation of riboflavin and its degradation products is easily achieved using Ultra Performance LC in less than four minutes, a vast improvement in speed and quality of data as compared to what could be previously achieved with traditional HPLC. Additionally, the use of ACQUITY UPLC photodiode array detection in series with single quadrupole mass spectrometry provides a powerful system platform for positive compound identification of known compounds. The use of UPLC/PDA/MS also aided in the initial identification of other related compounds in the mixture.

REFERENCES

1. J Chromatogr A. 2000 Jun 9;881(1-2):285-97.
2. J Chromatogr B Biomed Sci Appl. 2000 Feb 28;739(1):219-24.

Application NOTE

Waters

Sales Offices:

Austria 43 1 877 18 07

Australia 61 2 9933 1777

Belgium and Luxembourg
32 2 726 1000

Brazil 55 11 5543 7788

Canada 800 252 4752

China 86 10 8451 8918

Czech Republic 420 2 617 11384

Denmark 45 46 59 8080

Finland 358 9 506 4140

France 33 1 3048 7200

Germany 49 6196 400600

Hong Kong 852 29 64 1800

Hungary 36 1 350 5086

India 91 80 2837 1900

Ireland 353 1 448 1500

Italy 39 02 27 4211

Japan 81 3 3471 7191

Korea 82 2 820 2700

Mexico 52 55 5200 1860

The Netherlands 31 76 508 7200

Norway 47 6 384 6050

Poland 48 22 833 4400

Puerto Rico 787 747 8445

Russia/CIS 7 095 931 9193

Singapore 65 6278 7997

Spain 34 93 600 9300

Sweden 46 8 555 11 500

Switzerland 41 62 889 2030

Taiwan 886 2 2543 1898

UK 44 208 238 6100

US 1 800 252 4752

WATERS CORPORATION
34 Maple St.
Milford, MA 01757 U.S.A.
T: 508 478 2000
F: 508 872 1990
www.waters.com

Waters

For Complete Confidence

Waters, Empower, ZQ, UPLC, Ultra Performance LC, ACQUITY UPLC, and UPLC are trademarks of Waters Corporation.
All other trademarks are the property of their respective owners.
©2006 Waters Corporation Produced in the U.S.A. Jan. 2006 720001447EN KJ-PDF

