

SIMULTANEOUS ANALYSIS OF GHB AND ITS PRECURSORS IN URINE USING LC/MS/MS

Michelle Wood¹, Marleen Laloup², Nele Samyn², Michael Morris¹, Peter Batjoens³ and Gert De Boeck²

¹ Waters Corp., Manchester, UK; ²National Institute of Criminalistics and Criminology (N.I.C.C), Brussels, Belgium;
³Waters Corp., Brussels, Belgium.



Figure 1: GHB.

INTRODUCTION

Gamma-hydroxybutyrate (GHB) is a metabolite of gamma-aminobutyric acid (GABA) and plays the role of a central neurotransmitter and neuromodulator. Since GHB is a normal component of mammalian metabolism, it is present in all tissues of the body. Typical urinary GHB concentrations are < 10 mg/L^{1,2}. In some countries GHB is used clinically as an intravenous anaesthetic and as a treatment for narcolepsy, alcoholism and opiate withdrawal. Over the last few years, GHB has been gaining popularity amongst club-goers as the recreational drug (Figure 1) where it is taken for its ability to produce feelings of euphoria and to enhance sexuality³⁻⁵. As a result of its potent prosexual effects, GHB has also been increasingly implicated in drug-facilitated sexual assault^{6,7}. Ingestion of the chemical precursors of GHB i.e. gamma-butyrolactone (GBL) and 1,4-butanediol (1,4-BD) also results in similar physiological effects since they are rapidly converted to GHB in the body⁸. Raised awareness of the effects of these drugs and their potential for misuse, in addition to their ease of availability, has resulted in a dramatic increase in the demand for their analytical determination in both biological specimens and putative drug preparations. The purpose of this study was to develop and validate a sensitive LC/MS/MS procedure that would enable the simultaneous quantification of GHB, GBL and 1,4-BD in urine.

METHODS AND INSTRUMENTATION

Samples

Calibrators and quality control (QC) samples.

Control urine was spiked with GHB, GBL and 1,4-BD to yield a series of calibrators at the following concentrations; 0, 1, 2, 5, 10, 20, 50 and 80 mg/L. Low and high QC samples were prepared by spiking control urine with the drugs to yield concentrations of 4 and 40 mg/L, respectively.

Authentic Samples

One hundred and eighty two authentic human urine samples were collected from club-goers attending a post dance-club 'chill-out' venue and were the result of 2 separate raids by the Belgian Police Department. The samples were analysed for GHB and the precursors using the newly-developed LC/MS/MS procedure. For comparative purposes, the samples were also analysed for GHB using a routinely used GC/MS procedure.

Sample Preparation

Urine samples were diluted (1:20) with an internal standard solution (GHB-d6 and GBL-d6, at a concentration of 2 mg/L in deionised water).

LC Conditions

LC system:	Alliance® 2795
Column:	Atlantis® dC ₁₈ column (3 x 100 mm, 5 µm) at 35 °C
Mobile phases:	(A): 0.1% aqueous formic acid (B): methanol
Isocratic elution:	90:10 (A:B)
Flow rate:	200 µL/min.
Inj. volume:	20 µL

Mass Spectrometry Conditions

Mass spectrometer: Quattro microTM mass spectrometer
 Ionisation mode: ES⁺
 Capillary voltage: 3.5 kV
 Source Temperature: 120 °C
 Desolvation gas: Nitrogen at 700 L/Hr, 350 °C
 MS/MS: Collision gas (argon) at
 5×10^{-3} mbar

RESULTS AND DISCUSSION

Multiple reaction monitoring (MRM) transitions were determined for GHB, the precursors and the internal standards i.e. GHB-d6 and GBL-d6 (Table 1). Figure 2 shows some examples of product ion spectra.

A series of urine calibrators was prepared. Following preparation i.e. simple dilution, the samples were analysed using LC/MS/MS. Figure 3 shows the MRM chromatograms obtained following the analysis of a control urine sample and the same sample enriched with GHB, GBL and 1,4-BD.

Quantification was achieved by integration of the area under the specific MRM chromatogram. For GHB and GBL, responses were calculated in reference to the integrated area of their respective deuterated internal standards. For 1,4-BD the response was calculated by reference to that of GHB-d6. Linear responses were obtained for GHB and 1,4-BD over the range investigated (1-80 mg/L). GBL produced a linear response over the range 1-50 mg/L.

Precision, intra-assay and interassay variation (as % CV) were all found to be highly satisfactory at < 7%. Analytical recoveries ranged from 90-107% (Table 2).

Compound	Precursor ion (m/z)	Product ion (m/z)	Cone Voltage (V)	Collision energy (eV)
GHB	105	87	10	7
GHB-d6	111	93	10	7
GBL	87	45	25	15
GBL-d6	93	49	25	15
1,4-BD	91	73	12	5

Table 1: MRM transitions and conditions for the measurement of GHB, GBL, 1,4-BD and their deuterated internal standards.

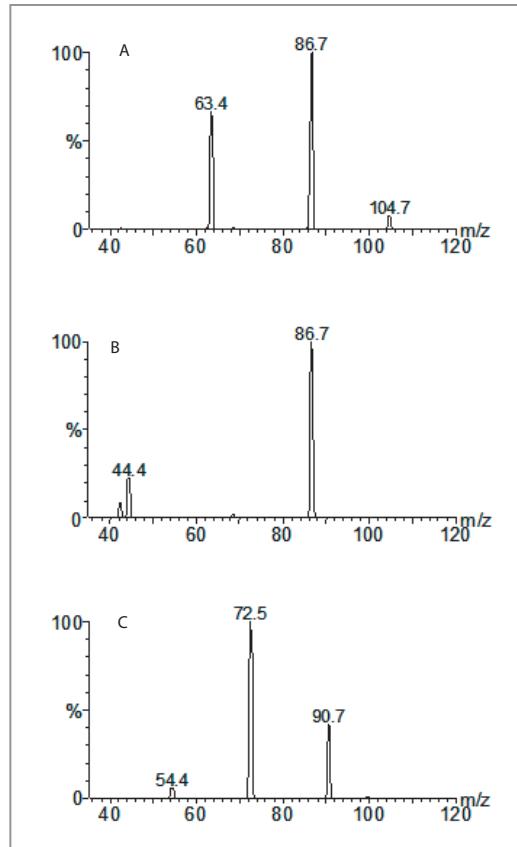


Figure 2: Product ion spectra for GHB (A), GBL (B) and 1,4-BD (C). Pure standards (5 mg/L) were infused into the mass spectrometer and the cone voltage (CV) optimised for the precursor ion*. CID was then performed and product ion spectra acquired under optimum conditions for the most abundant product ion.

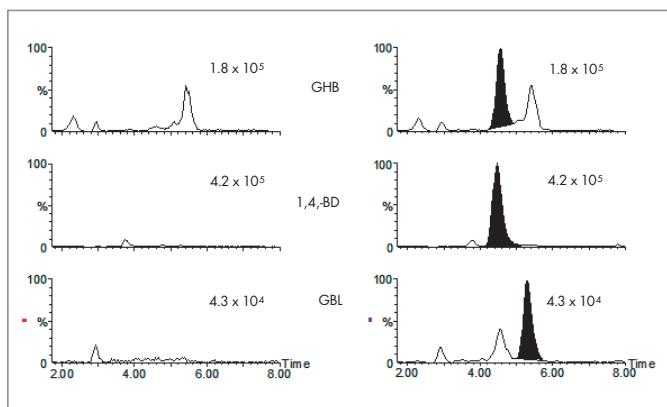


Figure 3: MRM chromatograms obtained with a single injection of a control urine sample (left-hand column) prepared by the dilution method and the same sample enriched with 10 mg/L of GHB, GBL and 1,4-BD (right-hand column). Peak intensity is shown in the top right-hand corner of each trace.

Compound	Low Control (4 mg/L)			High Control (40 mg/L)		
	Mean (mg/L)	Recovery (%)	%CV	Mean (mg/L)	Recovery (%)	%CV
Precision (n=5)						
GHB	4.0	100	3.0	41.0	103	0.5
GBL	3.8	95	4.2	40.2	101	3.9
1,4-BD	3.7	93	2.9	41.3	104	0.7
Intra-assay (n=5)						
GHB	3.9	98	3.2	42.7	107	3.5
GBL	3.7	93	3.2	36.1	90	2.9
1,4-BD	4.0	100	2.2	40.0	100	3.1
Intrarassay (n=5)						
GHB	4.1	103	5.3	40.0	100	3.4
GBL	4.0	100	6.6	39.8	100	6.3
1,4-BD	3.9	98	3.8	40.5	101	4.7

Table 2: Precision and analytical recovery data for GHB and its precursors in urine.

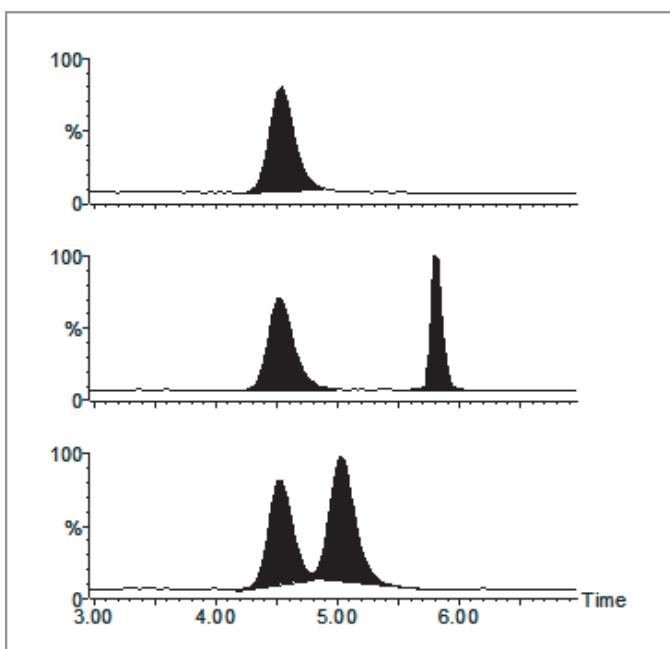


Figure 4: LC/MS analysis of the hydroxybutyric acid isomers. Ion chromatograms obtained following the analysis of gamma-hydroxybutyric acid (GHB) only (A) and GHB in the presence of alpha and beta-hydroxybutyric acid (traces B and C respectively). Peak intensity is shown in the top right-hand corner of each trace and is the sum of the responses obtained for both the protonated and the sodiated species i.e. m/z 105 + 127.

The limit of quantification was defined as the concentration of the lowest calibrator which was calculated to be with in $\pm 20\%$ of the nominal value and with a % CV less than 20%. For all of the analytes of interest, this criteria was met by the 1 mg/L calibrator and was sufficient to determine endogenous levels of GHB in the urine.

To investigate any potential interference in GHB quantification by its naturally occurring isomers i.e. alpha and beta-hydroxybutyric acid, standards were analysed using the developed LC/MS/MS method. Both compounds were shown to be chromatographically resolved from GHB and thus would not interfere in the quantification of the latter (Figure 4).

The utility of the LC/MS/MS method was demonstrated by the analysis of one hundred and eighty-two authentic urine samples. Seven samples contained GHB at concentrations > 2 mg/L. The same seven samples were independently identified by the more time-consuming, labour-intensive GC/MS method.

Only two, of these seven samples, were above the recommended interpretive cut-off concentration of 10 mg/L and were 956 mg/L and 1411 mg/L, respectively. These two samples were also positive for GBL. None of the authentic urine samples contained 1,4-BD.

CONCLUSIONS

The method presented here is the first demonstration of the use of LC/MS/MS for the simultaneous analysis of GHB and its precursors in urine samples. The method is simple and rapid (total analysis time of <12 min). It offers sufficient sensitivity to enable the measurement of endogenous levels of GHB and to identify exogenous ingestion.

The LC/MS/MS results obtained following the analysis of authentic samples, correlated with the more labour-intensive, time-consuming (~1 hour) GC/MS method.

The procedure offers several advantages over other published techniques;

1. It enables the simultaneous quantification of the GHB and the precursors in a single analysis; this can facilitate the identification of the chemical basis of any seized putative drug preparations or if present in the biological specimen, can provide information of the chemical nature of the ingested drug.
2. It involves fewer manipulations and is less time-consuming.

[APPLICATION NOTE]

Waters

THE SCIENCE OF WHAT'S POSSIBLE.TM

Although the data presented here indicate that the actual prevalence of GHB-positives might be quite low, the hype and publicity surrounding these drugs has led to a dramatic increase in the number of requests for their analysis in biological samples (and particularly in urine). The simplicity and speed of the described LC/MS/MS technique, should prove a useful means to meet this current increased demand on laboratories.

REFERENCES

1. Elliott SP. Gamma hydroxybutyric acid (GHB) concentrations in humans and factors affecting endogenous production. *Forensic Sci. Int* 2003;133:9-16.
2. LeBeau MA, Christenson RH, Levine B, Darwin WD and Huestis MA. Intra- and inter individual variations in urinary concentrations of endogenous gamma-hydroxybutyrate. *J. Anal. Toxicol* 2002;26:340-346.
3. Laborit H. Correlations between protein and serotonin synthesis during various activities of the central nervous system (slow and desynchronized sleep, learning and memory, sexual activity, morphine tolerance, aggressiveness and pharmacological action of sodium gamma-hydroxybutyrate). *Research Communications in Chemical Pathology and Pharmacology* 1972;3:51-81.
4. Ropero-Miller JD and Goldberger BA. Recreational drug current trends in the 90's. *Clin. Lab Med* 1998;18:727-746.
5. Bellis MA, Hughes K, Bennett A and Thomson R. The role of an international nightlife resort in the proliferation of recreational drugs. *Addiction* 2003;98:1713-1721.
6. ElSohly MA and Salamone SJ. Prevalence of drugs used in cases of alleged sexual assault. *J. Anal. Toxicol* 1999;23:141-146.
7. Ferrara SD, Frison G, Tedeschi L and LeBeau MA. Gamma-hydroxybutyrate (GHB) and related products. In: LeBeau MA and Mozayani A, eds. *Drug-Facilitated Sexual Assault (DFSA): A Forensic Handbook*. London: Academic Press, 2001:108-126.
8. Palatini P, Tedeschi L, Frison G, Padrini R, Zordan R and Orlando R et al. Dose-dependent absorption and elimination of Áhydroxybutyric acid in healthy volunteers. *Eur. J. Pharmacol* 1993;45:353-356.
9. Fieler EL, Coleman DE and Baselt RC. GHB concentrations in pre and post-mortem blood and urine [Letter]. *Clin. Chem* 1998;44:692.

Waters instruments are in vitro diagnostic devices compliant with the EU directive 98/79/EC and the United States Food and Drug Administration Quality System Regulations and current Good Manufacturing Practices. Waters Corporation is ISO 13485:2003 certified.

Austria and European Export (Central South Eastern Europe, CIS and Middle East) 43 1 877 18 07, Australia 61 2 9933 1777, Belgium 32 2 726 1000, Brazil 55 11 5094 3788, Canada 1 800 252 4752 x2205, China 86 10 8586 8899, CIS/Russia 7 095 336 7000, Czech Republic 420 2 617 1 1384 Denmark 45 46 59 8080, Finland 358 9 5659 6288, France 33 1 30 48 72 00, Germany 49 6196 400600, Hong Kong 852 29 64 1800 Hungary 36 1 350 5086, India and India Subcontinent 91 80 2837 1900, Ireland 353 1 448 1500, Italy 39 02 27 421 1, Japan 81 3 3471 7191 Korea 82 2 820 2700, Mexico 52 55 5200 1860, The Netherlands 31 76 508 7200, Norway 47 6 384 60 50, Poland 48 22 833 4400 Puerto Rico 1 787 747 8445, Singapore 65 6273 1221, Spain 34 93 600 9300, Sweden 46 8 555 11 500, Switzerland 41 62 889 2030 Taiwan 886 2 2543 1898, United Kingdom 44 208 238 6100

All other countries: Waters Corporation U.S.A. 1 508 478 2000/1 800 252 4752



Waters
THE SCIENCE OF WHAT'S POSSIBLE.TM

Waters, Atlantis, Quattro micro, The Science of What's Possible and Alliance are trademarks of Waters Corporation.
All other trademarks are acknowledged.
©2007 Waters Corporation. Printed in the U.S.A.
July 2007 720001548EN MC-PDF

Waters Corporation
34 Maple Street
Milford, MA 01757 U.S.A.
T: 1 508 478 2000
F: 1 508 872 1990
www.waters.com