Development of a Quantitative SPE/LC/MS/MS Assay for Multiple Amyloid Beta Peptides in Cerebrospinal Fluid

Erin Chambers¹, Mary E. Lame², Diane M. Diehl¹, Sarah Grimwood², and Yanhua Zhang³ ¹Waters Corporation, Milford, Massachusetts 01757, ²Pfizer Neuroscience Research Unit, Groton, CT 06340, ³Pfizer PDM, Groton, CT 06340

INTRODUCTION

IZEI

The deposition/formation of insoluble aggregates of amyloid β peptides (A β) in the brain is considered to be a critical event in Alzheimer's Disease (AD). Therapeutic strategies have focused on small molecule inhibitors or immunotherapy in lowering amyloid beta production or enhancing its clearance. Therefore, the need for highly sensitive and robust methods for quantitation of amyloid peptides in cerebrospinal fluid and their relationship to AD is of critical interest to many researchers. However, the analysis of these Aß peptides is extremely challenging due not only to the relatively low abundance in biological fluids but also because they have a propensity for aggregation, they are potentially bound by other proteins and have the tendency to form oligomers.

Measurement of these peptides routinely employs immunoassays (for their selectivity and sensitivity), or tedious immunoprecipitation followed by SPE. Immunoassays require more time to develop than LC/MS/MS assays, they require multiple assays for multiple AB peptides, and have a limited linear dynamic range relative to LC/MS/MS. Immunoassays are subject to cross reactivity and non-specific binding, require expensive antibodies and rely on the selectivity of the antibody to enrich the sample/species. They are labor intensive and assay imprecision and matrix interferences are common issues. Therefore, there exists a need for a high throughput selective bioanalytical method based on LC/MS/MS with sample preparation capable of recovering pg/mL levels of amyloid peptides in the presence of high concentrations of interfering proteins and peptides.

While the time required developing immunoassays may be acceptable in later stages of the drug development process, it is hardly practical for the earlier stages, where a high throughput, reliable method capable of quantitating multiple peptides in a class is desirable.

This work focuses on the development of LC, MS, and selective SPE sample preparation methods for the 1-38, 1-40, and 1-42 fragments of amyloid precursor protein (APP), in support of preclinical studies. The use of a single, high throughput assay for multiple Aß peptides, without time consuming immunoprecipitation steps, was developed and validated. The A β class of peptides, in particular, presents many unique analytical challenges including non-specific binding, poor solubility, aggregation, and lower MS sensitivity. Steps were taken at all stages of method development to minimize or eliminate the impact of these issues.

As strategies emerge for disease modification in AD, the guantification of multiple A β species, in addition to A β 38, 40 and 42, that may be linked to AD pathology may help to offer more insight to this disease and its progression. The method described herein shows promise for adaptation to quantitate those peptides as well.

Amyloid β 1-38, MW 4132 PI 5.2 DAEFRHDSGYEVHHQKLVFFAEDVGSNKGAIIGLMVGG **Amyloid β 1-40,** MW 4330 PI 5.2 DAEFRHDSGYEVHHQKLVFFAEDVGSNKGAIIGLMVGGVV **Amyloid β 1-42,** MW 4516 PI 5.2 DAEFRHDSGYEVHHQKLVFFAEDVGSNKGAIIGLMVGGVVIA

Figure 1: Amino acid sequence and pI data for amyloid β peptides 1-38, 1-40, and 1-42.

EXPERIMENTAL

	UPLC [®] Method Conditions
Column:	ACQUITY UPLC [®] BEH C ₁₈ 300Å 2.1 \times

Corannin	1.00				U 10	500/ CEI	- · ·	
	150	mm,	1.7 μr	n				
Mobile Pha	ases:	A: 0	.3% N	IH₄OH	(by	volume)) in	H_2O

Gradient:

B: 90% Acetonitrile, 10% mobile phase A hold 90% A for 1 min., ramp to 55% A in 5.5 min, hold 0.2 min, return to initial Flow Rate: 0.2 mL/min 10 µL

Injection: Temperature: 50 °C

MS Conditions

System: Waters Xevo[™] TO triple guadrupole mass spectrometer operated in ESI+ MRM mode

Desolvation Gas Flow: 800 L/hr

Source Temperature: 120 °C

Desolvation Temperature: 450 °C

Collision Cell Pressure: 2.6 X 10 (-3) mbar MRM Transitions and conditions:

see Table 1 Sample Pre-treatment

200 μ L CSF (human, monkey, or spiked artificial CSF + 5% rat plasma) was diluted 1:1 with 5M guanidine HCL and shaken at room temperature for 45 minutes. This was then diluted further with 200 μ L 4% H₃PO₄ in H₂O. Note: for spiked samples, samples were allowed to equilibrate at room temperature for 30 min. after spiking, prior to dilution with quanidine HCl.

Solid Phase Extraction (SPE) Oasis[®] MCX in µElution 96-well format

Condition: 200 µL methanol Equilibrate: 200 μ L 4% H₃PO₄ in H₂O Load: 600 µL pre-treated sample Wash 1: 200 µL 4% H₃PO₄ in H₂O Wash 2: 10% ACN in H_2O Elute: 2 X 25 µL 75:15:10 ACN:H₂O:conc. NH₄OH Dilute: 25 μ L H₂O Inject 10 µL

	Precursor	Product	Product	Cone	Collision
Peptide Name	lon	lon	lon ID	voltage (V)	energy (eV)
Amyloid β 1-38	1033.5	1000.3	b 36	33	23
Amyloid β 1-38 N15 IS	1046	1012.5		30	22
Amyloid β 1-40	1083	1053.6	b 39	33	25
Amyloid β 1-40 N15 IS	1096	1066.5		35	22
Amyloid β 1-42	1129	1078.5	b 40	28	30
Amyloid β 1-42 N15 IS	1142.5	1091.5		35	28

Table 1: MRM transitions and MS conditions for the amyloid β peptides and their N15 labeled internal standards



Figure 2: Representative ESI+ MSMS spectrum for amyloid β 1-42

TO DOWNLOAD A COPY OF THIS POSTER, VISIT WWW.WATERS.COM/POSTERS

RESULTS AND DISCUSSION

The greatest challenges encountered in the development of these methodologies were overcoming solubility, adsorption, and aggregation issues and obtaining adequate selectivity and sensitivity to meet the application requirements. Proper mobile phase and injection solvent composition, as well as judicious choice of SPE elution solvent are just a few of the factors that were key to addressing these problems.

MASS SPECTROMETRY

MS was performed in positive ion mode as CID of the 4+ precursor vielded several distinct product ions (representative spectrum shown in Figure 2) corresponding to inherently specific b sequence ions. MSMS in negative ion mode vielded a dominant water loss. Figure 3 demonstrates one example of the specificity difference between both methods. Although overall sensitivity was higher in solvent standards using the negative ion method, the sensitivity difference was mitigated in the presence of matrix and the improved specificity and signal to noise in positive ion mode proved critical for accurate quantitation in CSF samples.

UPLC

Separation of the three amyloid β peptides is shown in Figure 4. While exact % NH4OH in the mobile phase was critical for negative ion sensitivity, the signal in ESI+ proved to be more robust to subtle changes in mobile phase composition, providing a minimum of >24 hour LC/autosampler stability. In contrast, 50% or more of the ESI- signal was lost after 10-12 hours due to the natural change in NH_4OH concentration (volatility) in the mobile phase. This further reinforced the robustness of an ESI+ MS method.

SOLID PHASE EXTRACTION (SPE)

SPE was performed using Oasis[®] MCX, a mixed-mode sorbent, to enhance selectivity of the extraction. The sorbent relies on both reversed-phase and ion exchange retention mechanisms to selectively separate the amyloid β fraction from other high abundance polypeptides in complex CSF samples. The specific 96-well format used, Oasis[®] µElution, provided significant concentration while eliminating evaporation and reconstitution, minimizing peptide losses. In addition, binding of the peptides by ion-exchange imparted orthongonality into the overall method.

During initial method development a high degree of non-specific binding (NSB) was observed when artificial CSF was extracted. 5% rat plasma (having a different amyoid β sequence) was added to eliminate the NSB.

The SPE was one of the more critical aspects of the overall methodology. Very selective isolation of the amyloid fraction coupled to the resolution of standard flow UPLC enabled very rapid analysis of pre-clinical samples.

LINEARITY, ACCURACY AND PRECISION

N15 labeled internal standards were used for each peptide. Standard curves for all 3 amyloid β peptides were linear, with 1/xweighting from 0.1 to 10 ng/mL in artificial CSF + 5% rat plasma. A representative standard curve for amyloid β 1-38 is shown in Figure 5. Basal levels of the amyloid peptides were quantitated using both standard curves prepared from over-spiked human CSF and from artificial CSF + 5% rat plasma, and calculated basal levels were not statistically different. The artificial CSF was chosen as it is a less expensive and more readily available matrix. Basal levels of amyloid B 1-42 extracted from 3 sources of human and 1 source of monkey CSF are shown in Figure 6. Statistics for the determination of basal levels of all 3 amyloid β peptides are shown in Table 2.

Overspike QC samples were prepared in 3 sources of pooled human CSF and 1 source of pooled monkey CSF at 0.2, 0.8, 2, and 6 ng/mL. Accuracy and precision values met the regulatory criteria for LC/MS/MS assays. Representative results from QC sample analysis are shown in Table 3.



Figure 3: Comparison of MS specificity using either negative ion (water loss product ion) or positive ion (b sequence ion product ion) electrospray





Figure 4: UPLC/MS/MS analysis of amyloid β 1-38, 1-40 and 1-42 peptides extracted from artificial CSF + 5% rat plasma









Waters THE SCIENCE OF WHAT'S POSSIBLE.™

Figure 6: Representative chromatogram showing basal levels of amyloid β1-42 in 3 sources of human and 1 source of mon-

leta 1	-38				Amyloid Beta 1	-42				6
	Human CSF		Human CSF			Human CSF		Human CSF		
	Pool 1	Human CSF	Pool 3	Cyno CSF		Pool 1	Human CSF	Pool 3	Cyno CSF	
e #	ng/mL	Pool 2 ng/mL	ng/mL	Pool 1 ng/mL	Replicate #	ng/mL	Pool 2 ng/mL	ng/mL	Pool 1 ng/mL	
	1.585	2.354	1.014	1.713	1	0.519	0.616	0.421	0.675	
	1.650	2.103	1.371	1.605	2	0.421	0.656	0.481	0.621	
	1.614	2.464	0.950	1.947	3	0.542	0.644	0.534	0.623	
	1.657	1.939	1.608	1.541	4	0.471	0.567	0.348	0.659	
	1.820	2.158	1.471	1.675	5	0.476	0.573	0.487	0.700	_
	1.486	1.995	1.167	1.644	6	0.561	0.713	0.510	0.688	/
۱	1.635	2.169	1.264	1.688	Mean	0.498	0.628	0.463	0.661	
ation	0.110	0.204	0.262	0.140	Std. Deviation	0.052	0.055	0.068	0.033	
/	6.7	9.4	20.7	83	% C V	10.4	8.7	14.7	5.1	
leta 1	-40									

	Human CSE		Human CSE	
	Pool 1	Human CSF	Pool 3	Cvno CSF
e#	ng/mL	Pool 2 ng/mL	ng/mL	Pool 1 ng/ml
	3.083	4.031	2.541	3.699
	3.391	3.776	2.593	3.989
	3.292	3.598	2.580	3.525
	2.884	3.533	2.612	3.956
	3.131	3.230	2.508	3.284
	3.656	3.619	2.490	3.595
-	3.240	3.631	2.554	3.675
ation	0.269	0.266	0.049	0.268
/	83	73	19	73

Table 2: Baseline levels of amyloid β peptides in 3 sources of pooled human CSF and 1 source of pooled monkey CSF.

γloid β ´	1-38					
)ver-						
pike		Mean			Number of	
onc.		Calculated	Std.		Replicates	Mean
g/mL	QC Conc.	Conc.	Deviation	% CV	Passed	Accuracy
0.2	1.49	1.355	0.071	5	3/3	90.9
0.8	2.09	1.843	0.118	6	3/3	88.2
2	3.29	3.287	0.319	10	3/3	99.9
6	7.29	7.701	0.478	6	3/3	105.6

γloid	β	1-40
lvor		

spike		Mean			Number of	
Conc.		Calculated	Std.		Replicates	Mean
ng/mL	QC Conc.	Conc.	Deviation	%CV	Passed	Accuracy
0.2	2.75	2.359	0.015	1	2/3	85.8
0.8	3.35	3.054	0.016	1	2/3	91.2
2	4.55	3.929	0.011	0	3/3	86.4
6	8.55	8 209	0.5	6	3/3	96

Amγloid β 1-42

)ver- spike Sonc. g/ml	OC Conc	Mean Calculated	Std.	% CV	Number of Replicates	Mean
g/inc	ac conc.	CONC.	Deviation	70 C V	Fasseu	Accuracy
0.2	0.663	0.655	0.07	11	3/3	98.8
0.8	1.263	1.145	0.058	5	3/3	90.7
2	2.463	2.403	0.121	5	3/3	97.5
6	6.463	5.986	0.701	12	3/3	92.6

Table 3: Representative results from analysis of QC samples prepared in pooled human CSF, source 3

CONCLUSIONS

- 1. An SPE-LC/MS/MS bioanalytical method was developed and validated for the simultaneous quantitation of multiple amyloid β peptides in human and monkey CSF.
- 2. The combination of a highly selective extraction method based on mixed-mode SPE in µElution format and the resolution of UPLC chromatography was key to the accurate, precise and reliable quantitation of 3 major amyloid β peptides in human and monkey CSF^{\cdot}
- 3. The use of positive ion MSMS and b ion sequence fragments provided the MS specificity required for this application.
- 4.96 samples can be extracted and ready for injection in < 30 minutes, providing the sample prep throughput required for pre-clinical studies.
- 5. The method described herein eliminates timeconsuming immunoassays or immunoprecipitation steps for pre-clinical work.
- 6. The mass range and sensitivity of the Xevo TQ MS enabled the selection of higher m/z precursors for fragmentation and the choice of highly specific b ion fragments, resulting in increased signal to noise and overall improved specificity for the assay.
- 7. This approach also allows one assay for the measurement of several different amyloid β peptides simultaneously in one sample with selectivity, specificity and in a high-thoughput format while still achieving the high sensitivity required for low level endogeneous amyloid β peptides. This represents a significant benefit as an ELISA assay would require multiple assays with multiple antibodies.

SELECTED REFERENCES

- 1. T.A. Lanz, J.B. Schachter, Journal of Neuroscience Methods 169 (2008) 16-22.
- 2. T. Oe et al, Rapid Communications in Mass Spectrometry 20 (2006) 3723-3735.
- 3. JR Slemmon et al, Journal of Chromatography B 846 (2007) 24-31.
- 4. NT Ditto et al, Journal of Neuroscience Methods 182 (2009) 260-265.
- 5. T.A. Lanz, J.B. Schachter, Journal of Neuroscience Methods 157 (2006) 71-81.
- 6. MJ Ford at al, Journal of Neuroscience Methods 168 (2008) 465-474.
- 7. E. Portelius et al, Journal of Proteome Research, 6 (2007) 4433-4439.

ACKNOWLEDGEMENT

The authors wish to acknowledge Wenlin Li (PDM, Pfizer) for her early work on amyloid β peptides using immunoaffinity LC/MS/MS.

Waters Corporation 34 Maple Street, Milford, MA 01757 tel. (508) 478-2000, fax (508) 478-1990

http://www.waters.com