

Identifying Leachables and Extractables from Packaging Materials

Baiba Čabovska,¹ Douglas M. Stevens,¹ A. John Cunningham,² Arthur E. Bailey²

¹Waters Corporation, Milford, MA, USA

²Mannkind Corporation, Danbury, CT, USA

APPLICATION BENEFITS

- Facilitates the daunting task of identifying unknown compounds in any field that deals with structural elucidation, such as Pharmaceutical, Chemical Material, and Food industries.
- Provides a workflow for the systematic identification of extractables.
- The same workflow applies to either GC or UPLC with QToF.

WATERS SOLUTIONS

Xevo® G2 QToF Mass Spectrometer

Atmospheric Pressure Gas Chromatography (APGC)

MassLynx™ Software

MS^E Technology

MassFragment™ Software

KEY WORDS

Extractables, leachables, resins, monomers and oligomers, plasticizers, stabilizers, fillers, coloring agents, antioxidants, antistatic agents, elemental composition

INTRODUCTION

The Pharmaceutical industry is required by the U.S. FDA to demonstrate that no toxic or harmful substances migrate from packaging materials into a drug during its expected product shelf life.¹⁻⁵ Similarly, in the Food and Cosmetics industries, there is significant interest in the investigation of packaging leachables present in their products. By definition, extractables are compounds that are extracted from packaging or device components under controlled extraction conditions. Leachables are compounds that migrate from the packaging into the product during its normal shelf life. In the ideal case, leachables are a subset of extractables. If a thorough and accurate identification – or at least compound class identification of all potential contaminants is not performed, it can lead to product recall, financial losses, and/or brand alienation for the company.⁶

The initial investigation, called a controlled extraction study, involves some type of solvent extraction, typically a reflux, microwave, or supercritical fluid extraction.⁷ The solvents chosen must cover a wide range of polarities to ensure that non-polar and polar analytes are extracted. The analytical techniques employed for analyzing extracts must be comprehensive to cover as many analytes as possible including GC-FID-MS (volatiles) and LC-UV-MS (non-volatiles).⁵

The challenge with the compounds observed in a controlled extraction study is their identification. Resin manufacturers rarely provide a complete list of all the additives in polymers used for packaging. The original ingredients can degrade or undergo chemical changes during the manufacturing process. Also, the resin manufacturer may not be aware of possible contaminants present within the compounds. Typical extractables include monomers and oligomers from incomplete polymerization reactions; plasticizers, stabilizers, fillers, coloring agents, antioxidants, and antistatic agents, as well as their degradants. Additionally, residues from detergents and mold release agents that can be present on the resin after the molding process.

EXPERIMENTAL

Sample preparation


Samples were prepared by microwave extraction. The samples of polypropylene and nylon (2 g) were extracted in 10 mL of isopropanol for 3 h at 70°C. After the extraction the supernatant was transferred to the GC vials.

MS CONDITIONS

MS System:	Xevo G2 QTof with 7890A GC
Column:	HP1-MS, 30 m x 0.32 mm, 1.0 µm film
Carrier gas:	He at 2 mL/min
Temp.:	35 °C for 5 min, 20 °C/min to 320 °C, hold 20.75 min
Injection port:	300 °C
Injection type:	1 µL splitless, 1 min purge
Makeup gas:	N ₂ at 500 mL/min
Scan range:	50 to 1,000 Da
Collision ramp for MS ^E :	15 to 25 eV
Data management:	MassLynx v. 4.1 Software

Many of the analytes obtained from single quadrupole GC/MS data can be identified using commercially available libraries, such as NIST. However, a difficulty arises for volatiles analysis when the compound of interest is not listed in the library, or when the sensitivity of a single quadrupole MS is not sufficient for a positive identification. Therefore, additional techniques, such as Atmospheric Pressure Gas Chromatography (APGC) and Quadrupole Time-of-Flight (QTof) described in this application note, are beneficial.⁸ Due to the absence of libraries for LC/MS data accurate mass data would vastly facilitate the non-volatile analysis. For both volatile and semi-volatile analysis performed here, MS^E data, acquisition on a quadrupole time of flight mass spectrometer, with commercially available structural elucidation tools proves to be valuable for identification of the unknown compounds.

Workflow

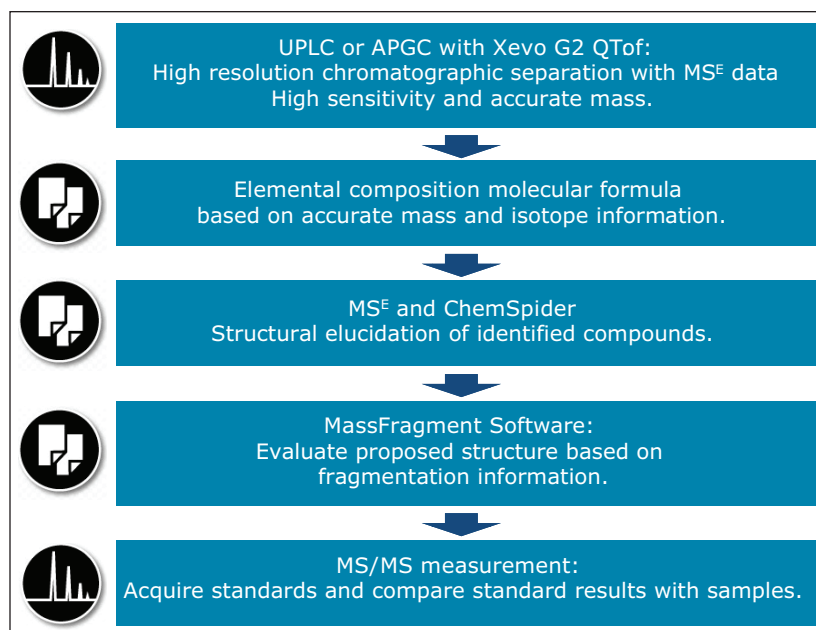

RESULTS AND DISCUSSION

Two widely available polymer materials were chosen for this study: polypropylene and nylon. In this application note, the identification of three different types of extractables is shown: an antioxidant, a monomer and a degradant of a monomer.

In the polypropylene sample, a peak (Peak A) was observed at a retention time of 26.3 min, as shown in Figure 1. Performing elemental composition analysis on the accurate mass APGC spectrum, shown in Figure 2, suggested a molecular formula of $C_{43}H_{63}O_3P$, as shown in Figure 3. The elemental composition software calculates the possible molecular formulas for the observed mass and also uses the isotope pattern algorithm to match the observed pattern with the theoretical one for each candidate molecular formula. In this case, there are two choices shown for the ion with the second being a closer match if only mass difference is considered. However, the combination of mass difference and isotope fit brings the correct one to the top of the list.

The APGC analysis was performed under dry source conditions,⁹ which promotes molecular ion (M^+) formation ahead of the protonated adduct ($[M+H]^+$). It is interesting to note that under high energy collision conditions the molecular ion fragments more easily than the protonated adduct; therefore the difference in the base peak was observed (646.4 versus 647.4) between the two channels, shown in Figure 2.

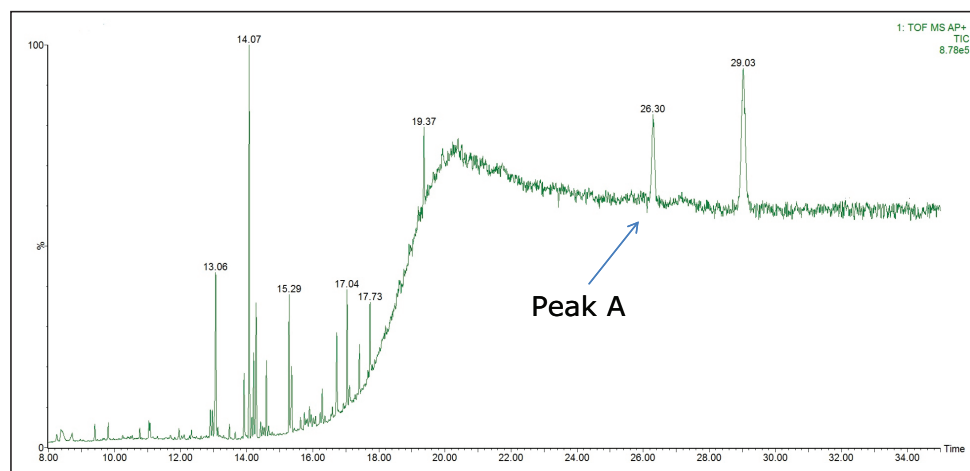


Figure 1. Polypropylene TIC.

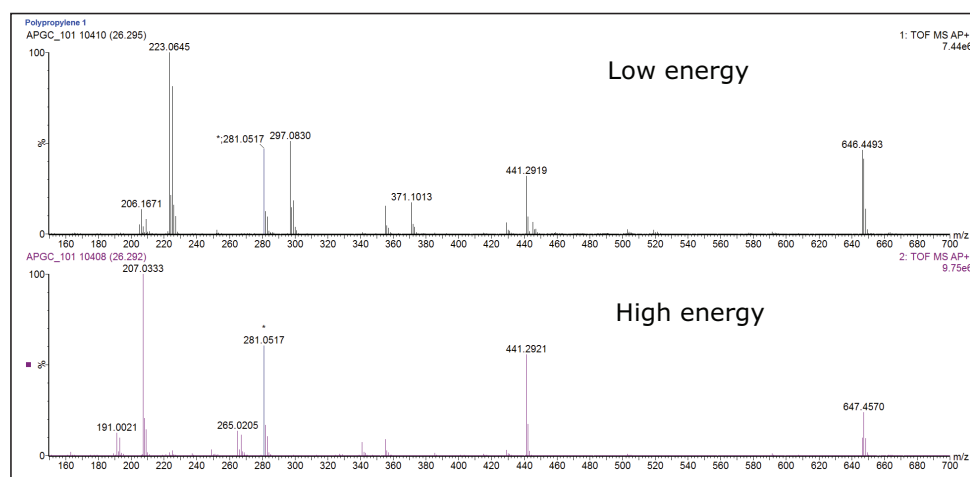


Figure 2. High and low energy spectra for Peak A.

Single Mass Analysis

Tolerance = 3.0 mDa / DBE: min = -1.5, max = 50.0

Element prediction: Off

Number of isotope peaks used for i-FIT = 3

Monoisotopic Mass, Odd and Even Electron Ions

265 formula(e) evaluated with 2 results within limits (up to 50 best isotopic matches for each mass)

Elements Used:

Mass	Calc. Mass	mDa	PPM	DBE	Formula	i-FIT	i-FIT Norm	Fit Conf %	C	H	O	P
646.4493	646.4515	-2.2	-3.4	12.0	C42 H63 O3 P	224.4	0.182	83.35	42	63	3	1
	646.4503	-1.0	-1.5	-1.0	C31 H66 O13	226.1	1.793	16.65	31	66	13	

Figure 3. Elemental composition data for Peak A.

Performing a search of the proposed elemental composition formula in ChemSpider gave Irgafos 168, shown in Figure 4, as the top answer when sorted by “# of References”, as described by Little, *et al.*¹⁰ Irgafos 168 is a trisarylphosphite processing stabilizer and protects the resin polymer, such as polypropylene, against oxidation during resin synthesis.

ChemSpider search results for $C_{42}H_{63}O_3P$. The search results are ordered by the number of references. The top result is Irgafos 168 (ID 82711), which has 38 references and 28 data sources. The second result is $C_{42}H_{63}O_3P$ (ID 2309060), which has 3 references and 3 data sources. The third result is $C_{42}H_{63}O_3P$ (ID 8028197), which has 2 references and 2 data sources.

ID	Structure	Molecular Formula	Molecular Weight	# of Data Sources	# of References
82711		$C_{42}H_{63}O_3P$	646.9216	28	38
2309060		$C_{42}H_{63}O_3P$	646.9216	3	3
8028197		$C_{42}H_{63}O_3P$	646.9216	2	2

Figure 4. ChemSpider search for $C_{42}H_{63}O_3P$, first match is Irgafos 168. The search hits are ordered by number of references and data sources.

Confidence in the identification was increased when another structural elucidation tool, Waters® MassFragment Software, was able to match several fragments observed in the high and low energy spectra to major fragment ions of Irgafos 168, as shown in Figure 5. MassFragment identifies bonds in precursor structure and then assigns a score based on the type and likelihood of the bond breakage. In addition, the number of bonds broken is listed. The lower the score (*e.g.* S:1.0, B:1.0 vs. S:4.5, B:2.0) the more probable the appearance of the fragment substructure.

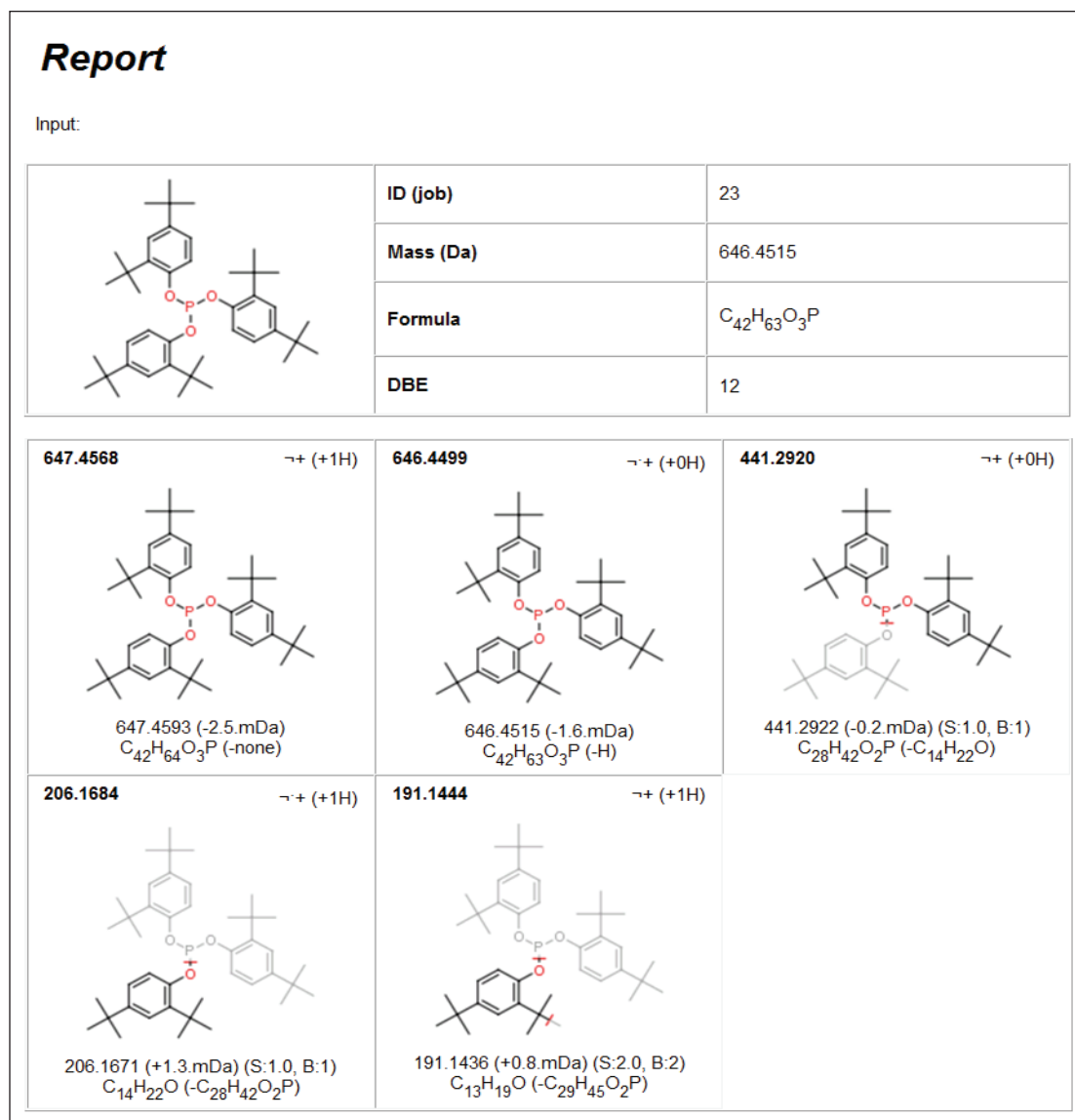


Figure 5. MassFragment Software report for confirmation of Irgafos 168.

The next step in this workflow is to purchase a standard and compare the retention time and fragmentation pattern with the sample.

Laurin lactam is a known starting material for the manufacturing of nylon. In the nylon extract the laurin lactam monomer (Peak B) is observed at a retention time of 15.93 minutes, as shown in Figure 6. The identity of the peak was confirmed by molecular formula and MassFragment following the workflow described in the previous example. A smaller peak is observed at a retention time of 16.07 minutes (Peak C). The measured mass is consistent with a molecular formula of $C_{12}H_{21}NO$, shown in Figure 7, which indicated that the peak was likely a laurin lactam degradant with an extra double bond in the molecule (laurin lactam monomer is $C_{12}H_{23}NO$). The parent ions in each spectra were confirmed by the presence of the in-source dimers ($2M+H$). For laurin lactam the observed dimer has m/z 395.3652 and for the degradant it is m/z 391.3324.

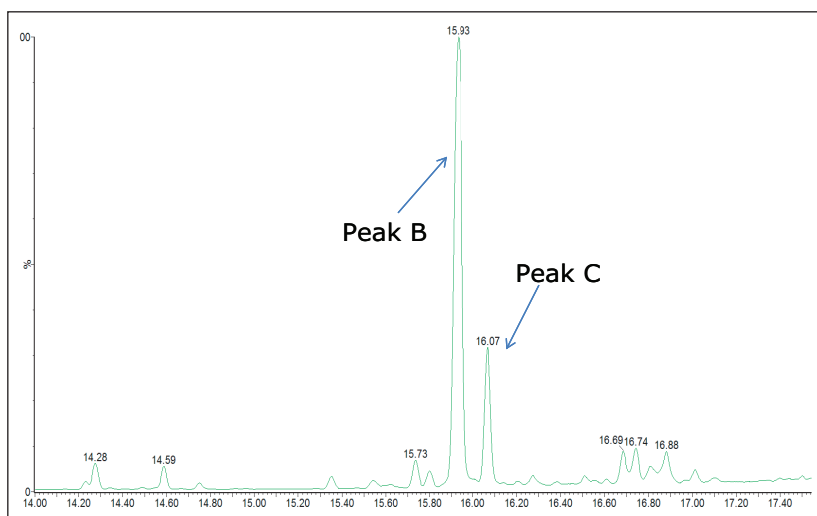


Figure 6. TIC for nylon extract.

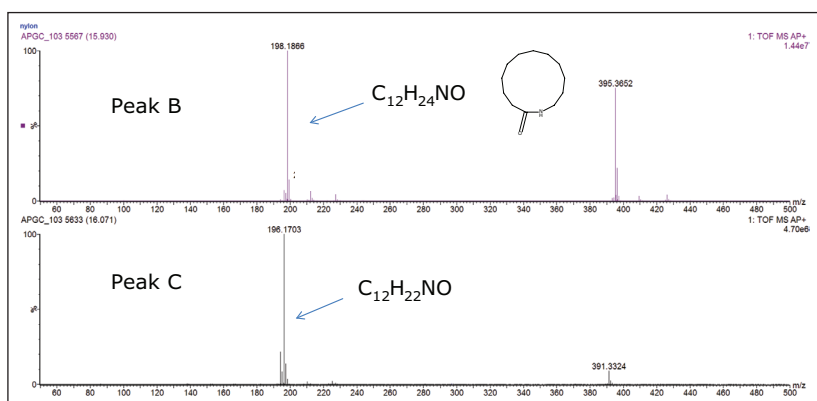


Figure 7. Spectra and molecular formula $[M+H]^+$ for Peaks B and C.

The ChemSpider search for $C_{12}H_{23}NO$ showed laurin lactam as the second top choice. The search of $C_{12}H_{21}NO$ did not provide any appropriate match based on the known compounds in the polymer.

Since a standard of this degradant is not likely to be available, the Xevo G2 QTof data allowed the assignment of a structure to this compound. It is not possible to determine the exact location of the double bond on the laurin lactam ring. However, in these types of studies it is not always necessary to determine an exact structure. It is sufficient if the compound's class has been identified. It was clear that the degradant is related to laurin lactam, therefore its toxicological profile was expected to be similar.

CONCLUSIONS

- Xevo G2 QTof is a valuable tool in the identification and structural elucidation of extractables. MS^E functionality allows simultaneous acquisition of precursor and fragment ions. Accurate mass and fragmentation information assists in the assignment of structures for many unknown compounds.
- Elemental composition and Mass Fragment Software provide the analyst with additional resources in cases when compounds of interest are not found in commercially available libraries.
- The workflow described can facilitate the daunting task of identifying the unknowns in any field that deals with structural elucidation, such as Pharmaceutical, Chemical Material, and Food industries.
- The fragments, the most likely molecular formula, and some chemical intuition based on ingredients known to be present can often provide a likely structure. In the extractable field a likely structure is often sufficient since the goal is to establish a safety threshold.

References

1. MP Balogh. "Testing the Critical Interface: Leachables and Extractables". LCGC, June 1, 2011.
2. F Mofatt. Extractables and Leachables in Pharma-Serious Issue. http://www.solvias.com/sites/default/files/solvias_whitepaper_web.pdf
3. Containers Closure Systems for Packaging Human Drugs and Biologics; Guidance for Industry. U.S. Department of Health and Human Services Food and Drug Administration Center for Drug Evaluation and Research (CDER) and Center for Biologics Evaluation and Research (CBER), Rockville, MD. May 1999.
4. D L Norwood, Q Fenge. Strategies for the analysis of pharmaceutical excipients and their trace level impurities. Am Pharm Rev. 7(5): 92, 94, 96-99, 2004.
5. Safety Thresholds and Best Practices for Extractables and Leachables in Orally Inhaled and Nasal Drug Products, Leachables and Extractables Working Group. Product Quality Research Institute (PQRI), 2006. (WWW.PQRI.ORG).
6. <http://www.webmd.com/pain-management/news/20110629/new-tylenol-recall-due-to-musty-odor>
7. M Ariasa, I Penichet, F Ysambert, R Bauza, M Zougaghc, Á Ríos. Fast supercritical fluid extraction of low- and high-density polyethylene additives: Comparison with conventional reflux and automatic Soxhlet extraction. J. of Supercritical Fluids, 50: 22-28, 2009.
8. Detection and Identification of Extractable Compounds from Polymers. Waters Technology Brief, no. 720004211en, January, 2012.
9. Determination of High Molecular Weight Phthalates in Sediments Using Atmospheric Pressure Chemical Ionization GC/MS. Waters poster, no. PSTRI34667160.
10. J L Little, A J Williams, A Pshenichnov, V Tkachenko. Identification of "known unknowns" utilizing accurate mass data and ChemSpider. JASMS. 23(1): 179-185, 2012.

Waters

THE SCIENCE OF WHAT'S POSSIBLE.™

Waters, Xevo, and UPLC are registered trademarks of Waters Corporation. MassLynx, MassFragment, and The Science of What's Possible are trademarks of Waters Corporation. All other trademarks are the property of their respective owners.

©2012 Waters Corporation. Produced in the U.S.A.
June 2012 720004391en AG-PDF

Waters Corporation
34 Maple Street
Milford, MA 01757 U.S.A.
T: 1 508 478 2000
F: 1 508 872 1990
www.waters.com

