

Increasing Efficiency of Method Validation for Metoclopramide HCl and Related Substances with Empower 3 MVM Software

Margaret Maziarz, Mark Wrona, and Sean M. McCarthy
Waters Corporation, Milford, MA, USA

APPLICATION BENEFITS

- Automated method validation workflow
- Reduced time to complete the steps required to test and document a validated method
- Compliance with regulations on data security, different user privileges, audit trails, data traceability, and electronic signature sign-off requirements

WATERS SOLUTIONS

ACQUITY UPLC® H-Class System

Empower® 3
Chromatography Data Software

Empower 3
Method Validation Manager (MVM)

ACQUITY UPLC Columns

ACQUITY® QDa® Detector

KEY WORDS

UPLC, method validation, metoclopramide HCl, sub-2-µm particles, mass detection, ApexTrack™ peak integration

INTRODUCTION

Method validation, which demonstrates that a method is suitable for its intended purpose,^{1,2} is an important regulatory requirement for pharmaceutical organizations and their supporting contract partners. A compliant laboratory must provide documented evidence and assurance that the analytical method used for testing a drug product's identity, quality, purity, and potency generates accurate and reliable results.

The validation process of an analytical method is a complex and demanding activity, consisting of many time-consuming steps. Some of these steps include creation of validation protocols, experimental work, reviewing and processing data, performing calculations, approving, and final reporting. Since some of these steps are prone to errors, a well-organized plan is essential for successful validation of an analytical method and to ensure that the appropriate regulations and guidelines are being followed.

Once validation is executed, adherence to the validation plan and specification is a critical compliance requirement. Any validation results not meeting the specifications must be clearly identified and addressed during the validation process. Raw and processed data must be appropriately stored and traceable by providing data security, audit trails, and automatic data documentation required for reviews and audits.

In this application note, we present validation of a UPLC® method for metoclopramide HCl and related substances using Empower 3 Method Validation Manager (MVM), an option for Empower 3 Chromatography Data Software.

We show how Empower 3 MVM tracked every step of the method validation process, identifying the steps and data that did not meet defined validation requirements.

Overall, we demonstrate that Empower 3 MVM automates the method validation workflow within a single software environment, reducing time and ensuring conformance to the validation requirements and acceptance criteria defined in the protocol.

EXPERIMENTAL

Method conditions

LC system: ACQUITY UPLC H-Class
 Column: ACQUITY UPLC CSH C₁₈,
 2.1 x 50 mm, 1.7 µm
 (Waters p/n 186005296)

Column temp.: 45 °C

Injection volume: 1.0 µL

Flow rate: 0.6 mL/min

Solvent A: 125 mM Formic acid
 in water

Solvent B: Water

Solvent C: Methanol

Separation: Gradient

Step	Time (minutes)	Solvent A (%)	Solvent B (%)	Solvent C (%)
1	Initial	10	85.0	5.0
2	5.0	10	30.0	60.0
3	5.5	10	30.0	60.0
4	5.6	10	85.0	5.0
5	7.5	10	85.0	5.0

Purge/Sample wash: 50:50 water/methanol

Seal wash: 90:10 water/acetonitrile

UV detector: ACQUITY UPLC PDA

UV setting: 210-400 nm
 (derived at 270 nm)

Mass detector: ACQUITY QDa
 (Extended Performance)

Ionization mode: ESI+, ESI-

Acquisition range: 100 – 440 *m/z*

Sampling rate: 10 pts/sec

Data: Centroid

System control,
 Data acquisition,
 and analysis: Empower 3 FR2
 CDS Software

RESULTS AND DISCUSSION

UPLC method for metoclopramide HCl and related substances

The UPLC method validated in this study was developed using a systematic method development protocol.³ An example of the UPLC chromatographic method for metoclopramide and related compounds is shown in Figure 1.

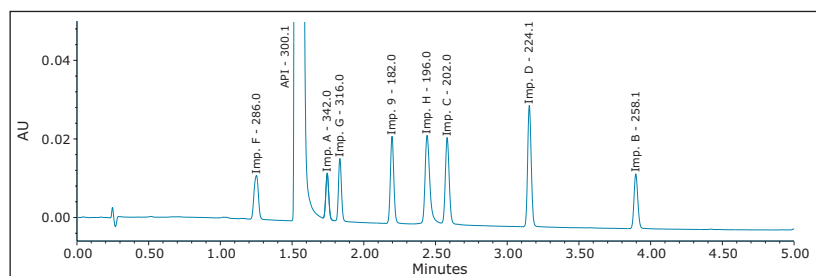


Figure 1. UPLC method for metoclopramide and related substances with UV at 270 nm.

About Empower 3 MVM

Empower 3 Method Validation Manager (MVM) is software that automates the validation process and enables efficient validation of chromatographic methods to ensure compliance to the validation requirements defined by the analytical laboratory. As shown in Figure 2, a validation workflow consists of many steps. A validation protocol (Figure 3) is created and used to execute the study. Once executed, Empower 3 MVM checks data for adherence with the validation requirements and flags any results that do not meet specifications. Validation results can be displayed in a report using validation report templates specific for each test available in Empower MVM software. The report templates can be customized as needed.

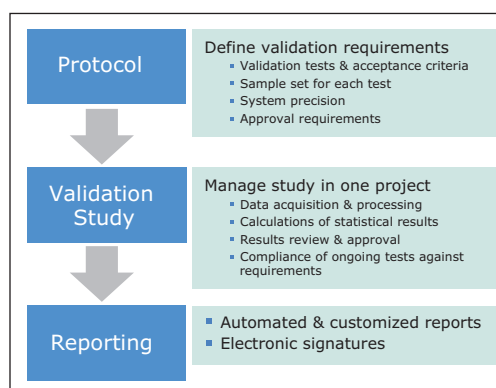


Figure 2. Workflow for validation of chromatographic methods with Empower 3 MVM.

Validation Protocol - Metoclopramide Related Sub in Validation_Metoclop_MVM_Demo as Lab_Manager/Admin...

File Edit View Tools Help

Method Classification

Compound Type: Drug Product

Analytical Method Type: Assay

Development Phase: Phase II

Protocol Comments: Protocol Template

Method/Study Approvals

Validation Protocol Approval: Jane Smith (Lab_Manager)

Validation Tests | System Precision | Approvals/Sign Offs

Validation Tests

Test Name	Test Description	Required	Test Status
1 Linearity	Linearity: 0.1 - 5.0 ug/mL	<input checked="" type="checkbox"/>	Sample Sets Incomplete
2 LDL/LQL Determination	Detection & Quantitation Limits	<input checked="" type="checkbox"/>	Sample Sets Incomplete
3 Accuracy	%Recovery - spiked drug tablet	<input checked="" type="checkbox"/>	Sample Sets Incomplete
4 Repeatability	6 preps at 0.1% level	<input checked="" type="checkbox"/>	Sample Sets Incomplete
5 Intermediate Precision	Inter Prec 1: Analysts 1 & 2	<input checked="" type="checkbox"/>	Sample Sets Incomplete
6 Specificity	Specificity	<input checked="" type="checkbox"/>	Sample Sets Incomplete
7 Robustness	Robust: temp, flow, wavelegth	<input checked="" type="checkbox"/>	Sample Sets Incomplete

Figure 3. Validation protocol method created within Empower 3 MVM project. The validation tests, acceptance criteria for each validation test, and requirements for approval are defined in the validation protocol method. The validation protocol is approved by a lab manager via electronic signature.

Method validation

We used an established validation protocol method approved by a lab manager (via electronic signatures) to execute the validation of our UPLC method for related substances of metoclopramide HCl in drug tablet formulation. The validation tests included linearity, detection and quantitation limits, accuracy, repeatability, intermediate precision, specificity, and robustness. System precision was evaluated for each validation test using five replicate injections of the sample, as recommended in the USP General Chapter <621> on Chromatography.⁴ The system precision criteria include:

- %RSD of retention times: $\leq 1.0\%$
- %RSD of peak areas: $\leq 2.0\%$
- USP resolution: ≥ 1.5
- Peak tailing: ≤ 1.5

1. Linearity

Method linearity for related substances was evaluated by analyzing seven concentrations of standard solutions ranging from 0.1 to 5.0 µg/mL. These concentrations corresponded to 0.02, 0.05, 0.1, 0.25, 0.5, 0.75, and 1.0% of the metoclopramide HCl target concentration of 0.5 mg/mL. We used Empower 3 MVM to calculate regression equation and correlation coefficients for a plot of average peak areas against the concentrations. Method linearity results generated by the software are displayed in Figure 4. The method shows linear relationship between the peak areas and concentrations for all related compounds with the correlation coefficients (r^2) greater than 0.999.

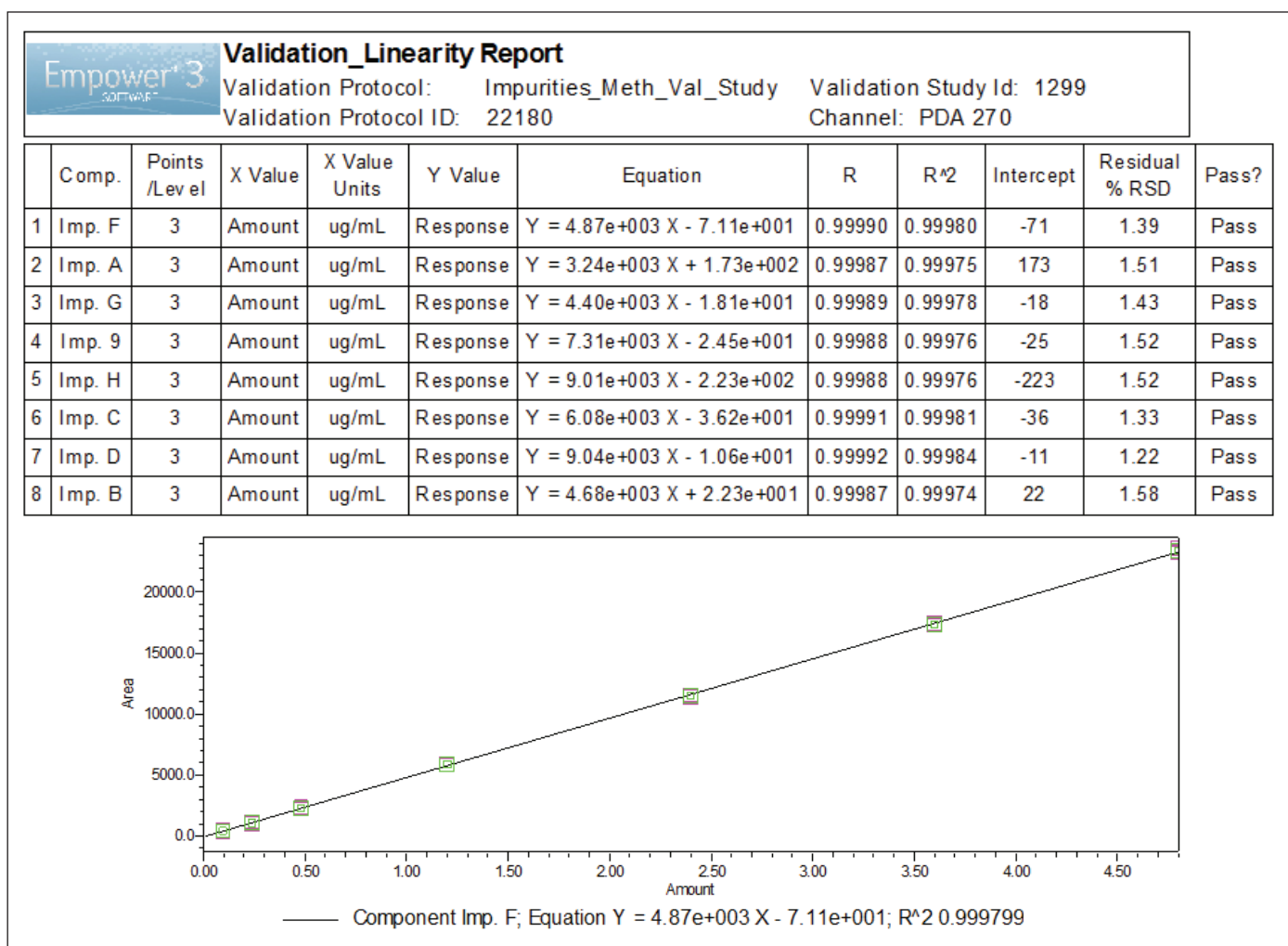


Figure 4. Method linearity results for metoclopramide related substances generated by Empower 3 MVM.

3. Accuracy

The accuracy of an analytical method includes quantitative determination of impurities in the presence of sample matrix components. Accuracy of our method was assessed by analyzing triplicate preparations of the drug tablet solutions spiked with related substances at 0.1, 0.5, and 1.0% levels in the presence of metoclopramide HCl concentration of 0.5 mg/mL. Accuracy results are summarized in Figure 7. The % recovery for all nine determinations ranged from 97 to 101% with %RSD \leq 4.21%, which passes the acceptance criteria of 90–110% and %RSD \leq 10%, respectively.

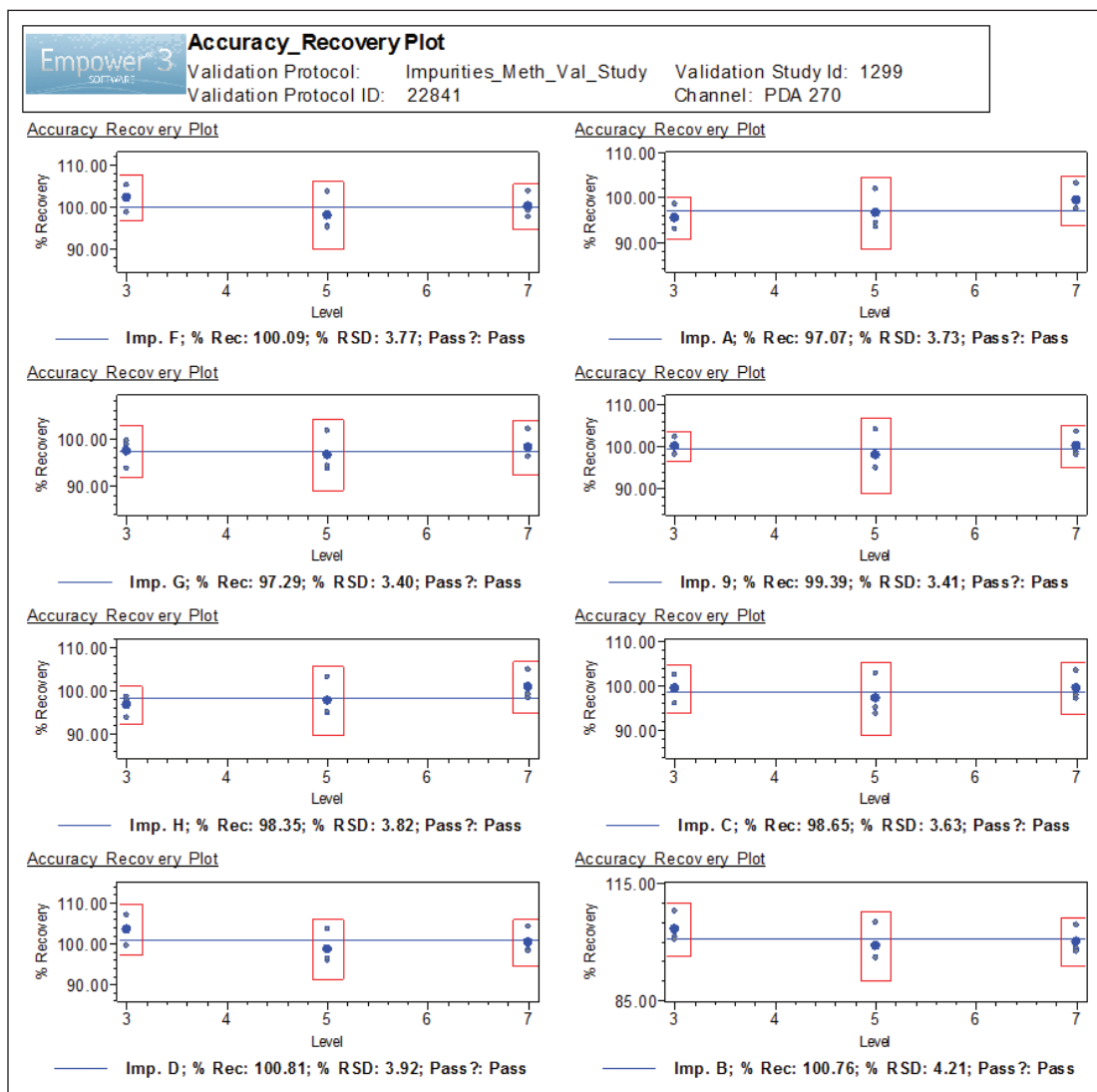


Figure 7. Accuracy results generated by Empower 3 MVM. Summary of 9 determination, 3 preparation at each levels: 0.1%, 0.5%, and 1.0%.

4. Repeatability

Method repeatability is a quantitative analysis of impurities from six independent preparations of the drug product by a single analyst. Repeatability of our method was demonstrated by spiking drug tablet sample solution with related substances at 0.1% level with respect to the metoclopramide HCl (API) concentration of 0.5 mg/mL. Repeatability results generated by analyst A (Figure 8) met the criteria for % recovery of $100 \pm 10\%$ and %RSD $\leq 10\%$.

5. Intermediate precision

Intermediate precision was evaluated by a different analyst, on a different day, using different instrument and column. Summary of results from six individual spiked drug tablet sample preparations generated by analysts A and B are shown in Figure 8. Overall, the intermediate precision results met the criteria for % recovery of $100 \pm 10\%$ and %RSD $\leq 10\%$.

<div> <div>Empower 3</div> <div> Intermediate Precision Validation Results Validation Protocol: Impurities_Meth_Val_Study Validation Study Id: 1299 Validation Protocol ID:11903 Channel: PDA Ch1 270nm@4.8nm, PDA </div> </div>						
Component: Imp. A Validation Result Id: 9208						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. A	1.792	Analyst A	6	94.8	2.31	2.44
2 Imp. A	1.825	Analyst B	6	96.6	2.05	2.13
Component: Imp. G Validation Result Id: 9209						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. G	1.890	Analyst A	6	97.3	3.45	3.54
2 Imp. G	1.922	Analyst B	6	96.8	1.60	1.66
Component: Imp. 9 Validation Result Id: 9210						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. 9	2.248	Analyst A	6	99.8	2.21	2.21
2 Imp. 9	2.305	Analyst B	6	98.8	1.17	1.18
Component: Imp. F Validation Result Id: 9285						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. F	1.311	Analyst A	6	101.4	2.62	2.58
2 Imp. F	1.343	Analyst B	6	101.8	1.40	1.38
Component: Imp. C Validation Result Id: 9212						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. C	2.636	Analyst A	6	99.5	2.05	2.06
2 Imp. C	2.696	Analyst B	6	96.9	0.89	0.92
Component: Imp. D Validation Result Id: 9213						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. D	3.186	Analyst A	6	102.4	3.28	3.20
2 Imp. D	3.254	Analyst B	6	97.4	1.15	1.18
Component: Imp. B Validation Result Id: 9214						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. B	3.914	Analyst A	6	102.2	3.35	3.28
2 Imp. B	3.996	Analyst B	6	100.0	2.21	2.22
Component: Imp. H Validation Result Id: 9211						
Component	RT Mean (min)	Experiment Group	Points	%Recovery Mean	Std. Dev.	%RSD of %Recovery
1 Imp. H	2.489	Analyst A	6	96.5	2.57	2.66
2 Imp. H	2.574	Analyst B	6	102.0	1.46	1.43

Figure 8. Repeatability (analyst A) and intermediate precision (analysts A and B) results generated by Empower 3 MVM.

6. Specificity

For the impurity test, specificity demonstrates that impurities can be separated and accurately measured in the presence of the sample matrix. This is typically done by spiking a drug substance or drug product with appropriate levels of impurities. In addition to demonstrating robust and reliable separation, it is important to identify that the desired components are not subject to interference with other species present in the sample. The UV peak purity determination is often used to show homogeneity of the chromatographic peak.

Specificity of our method was demonstrated by spiking drug tablet samples containing 0.5 mg/mL of metoclopramide HCl with related substances at 0.1% level. Accuracy and repeatability results show acceptable recoveries of each related substance. To demonstrate that the related substances are not coeluting with other components of the sample matrix, we used UV in conjunction with the MS spectral data as shown in Figure 9. Peak homogeneity was assessed using UV peak purity plot (Figure 9B). The peak purity angle is below the threshold angle, indicating the Impurity A peak is spectrally homogeneous. The mass spectral data provided additional information at the leading, apex, and tailing regions of the peak to confirm that only one mass is detected under the UV peak. The MS spectrum (Figure 9C) at the leading and tailing edge of the peak indicates the presence of an ion with mass of 342.0 m/z, which is specific to Impurity A. Overall, the UV peak purity plot and the MS spectrum shows that Impurity A is not coeluting with other peaks.

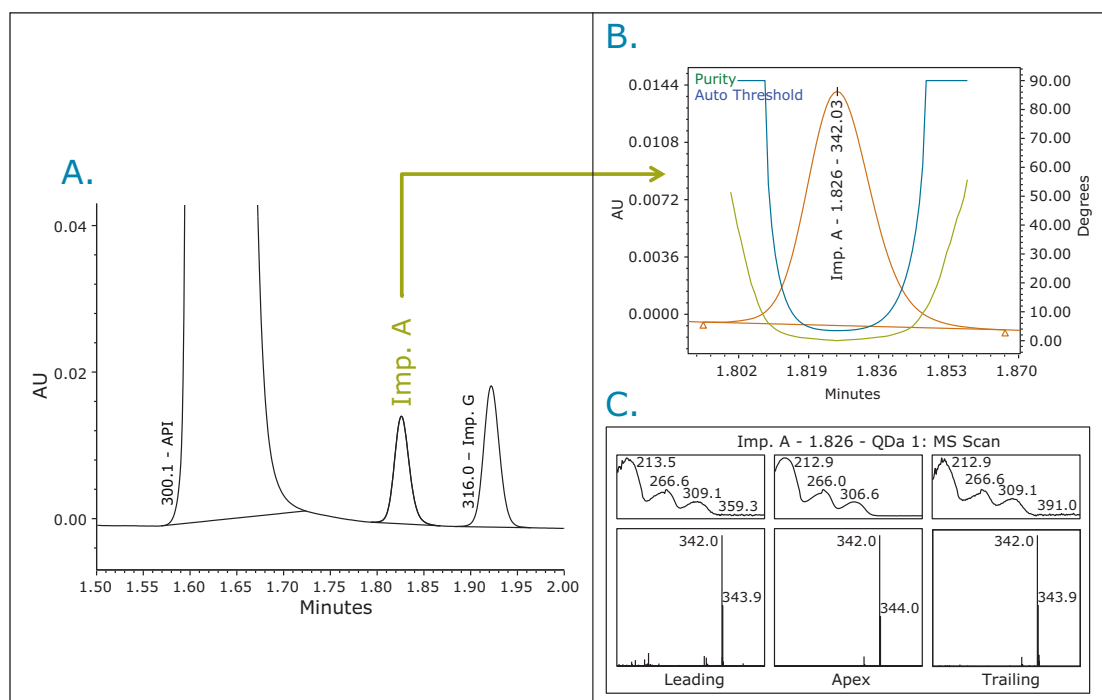


Figure 9. Peak homogeneity determination for specificity.

A. Accuracy sample with UV at 270 nm.

B. Peak purity plot of Impurity A.

C. UV and Mass profile of an Impurity A at the leading, apex, and tailing edge of the peak.

7. Robustness

Robustness is determined by the ability of the method to remain unaffected by the changes of chromatographic parameters. The parameters assessed in our study included:

- Column temperature: 45 ± 2.0 °C
- Flow rate: 0.6 ± 0.05 mL/min
- Wavelength: 270 ± 2 nm

Robustness was performed using a full factorial experimental design to investigate combination of different instrument conditions on the resolution between all the peaks, with a goal of achieving a minimum resolution of ≥ 2.0 for each peak. The robustness results in Figure 10 showed that the method met the criteria for resolution.

<div> <div>Empower 3 SOFTWARE</div> <div> Validation_Robustness_Report Validation Protocol: Impurities_Meth_Val_Study Validation Study Id: 1299 Validation Protocol ID: 22535 Channel: PDA Ch1 270nm@4.8nm </div> </div>									
Assessed Field: USP Resolution									
	Validation Result Id	Component	RT Mean (min)	Experiments	Assessed Field	Mean Rs	Lower Limit of Mean Rs	Upper Limit of Mean Rs	Pass /Fail
1	17451	API	1.617	8	USP Resolution	4.8	4.7	4.9	Pass
2	17487	Imp. A	1.825	8	USP Resolution	4.0	3.9	4.2	Pass
3	17489	Imp. G	1.923	8	USP Resolution	2.8	2.4	3.1	Pass
4	17491	Imp. 9	2.312	8	USP Resolution	10.3	10.1	10.5	Pass
5	17493	Imp. H	2.597	8	USP Resolution	6.2	5.5	7.0	Pass
6	17495	Imp. C	2.712	8	USP Resolution	2.4	2.0	2.8	Pass
7	17497	Imp. D	3.267	8	USP Resolution	12.9	12.3	13.4	Pass
8	17499	Imp. B	4.006	8	USP Resolution	17.1	16.6	17.6	Pass

Figure 10. Robustness results. Resolution for each component was ≥ 2.0 .

CONCLUSIONS

We successfully validated the UPLC method for related substances of metoclopramide HCl for linearity, detection and quantitation limits, accuracy, repeatability, intermediate precision, specificity, and robustness using Empower 3 MVM. Validation results showed that the method is linear, accurate, repeatable, precise, specific, and robust for all related substances tested in this study. The quantitation limit of all related compounds was below the reporting threshold of 0.1% or 0.5 µg/mL.

Empower 3 MVM software streamlined the entire validation process, from creating a validation protocol method to acquiring, reviewing, analyzing, approving, and reporting validation data. The updated tool/message center provided status of each validation test in a study and whether the results met the acceptance criteria, and flagged any out-of-specification results. Using ApexTrack for automated peak integration enabled consistent integration of all peaks during the validation process. Using the ACQUITY QDa Detector for mass detection in conjunction with UV detection enabled quick determination of peak purity using mass and UV

spectral data. Finally, the validation results and validation study were reviewed and approved via electronic signatures.

Empower 3 MVM is compliant-ready software that can be easily adapted by any analytical laboratory to improve efficiency of the chromatographic method validation process and to ensure conformance to validation requirements.

References

1. ICH Q2B, Validation of Analytical Procedures, International Conference on Harmonization, November 1996.
2. USP General Chapter, <1225>, Validation of Compendial Procedures, The United States Pharmacopeia Convention, official May 1, 2013.
3. Maziarz M, McCarthy SM, Wrona M. Improving Effectiveness of Method Development for Metoclopramide HCl and Related Substances Using a Systematic Screening Protocol. Waters Application Note; 2014: 720005026en.
4. USP General Chapter, <621>, Chromatography, USP36-NF31, The United States Pharmacopeia Convention, official December 2013.

Waters

THE SCIENCE OF WHAT'S POSSIBLE.®

Waters, The Science of What's Possible, ACQUITY, ACQUITY UPLC, UPLC, QDa, and Empower are registered trademarks of Waters Corporation. ApexTrack is a trademark of Waters Corporation. All other trademarks are the property of their respective owners.

©2014 Waters Corporation. Produced in the U.S.A. July 2014 720005111EN AG-PDF

Waters Corporation
34 Maple Street
Milford, MA 01757 U.S.A.
T: 1 508 478 2000
F: 1 508 872 1990
www.waters.com