

# Using LC/MS/MS 6410 for Analysis of Chloramphenicol, Thiamphenicol, and Florfenicol in Fish Samples

# **Application Brief**

Xiaorong Ran, Tao Bo

Chloramphenicol is a banned compound in the EU. It is a zero tolerance compound, and some methods have already been developed for this analysis. The LC/MS/MS is often used for its greater sensitivity and higher selectivity.

In many cases, not only chloramphenicol but also thiamphenicol and florfenicol are also found. This method uses a simple method for detecting all three compounds within less than 6 minutes. Furthermore, the results gave the good results, showing the performance at the negative mode.

# **Experimental**

#### **LC Conditions**

Column Agilent ZORBAX Eclipse Plus, 2.1 mm  $\times$  50 mm, 1.8  $\mu$ m

Mobile phase A: Water

B: Methanol

Flow rate 0.4 mL/min

Gradient 0–2 min/B 10% to 90%; 2–3 min/B 90%; 3.01/B 10%

 $\begin{array}{lll} \text{Stop time} & \text{6 min} \\ \text{Column compartment temperature} & \text{45 °C} \\ \text{Injection} & \text{5 } \mu \text{L} \\ \end{array}$ 

MSD Condition

Instrument Agilent 6410A triple quadrupole LC/MS system

Source ESI –

# **Highlights**

- Using a simple RRLC method can separate the compounds well within 6 minutes
- Quite high sensitivity in negative mode
- The ISTD method can remove the matrix effect and minimumize the sample preparation interference



#### **Sample Preparation**

All SPE cartridges are conditioned with 2 mL of water before use.

- 1. Honey, 1 g sample is diluted to 5 mL with water and 25 µL 10 ppb IS is added. The solution is loaded onto the SPE cartridge and allowed to stand for 5 min. Elution is performed with 10 mL ethyl acetate. The eluate is collected and the solvent is evaporated under a nitrogen stream at 40 °C. The residue is redissolved in 1 mL methanol and put in an ultrasonic bath for 1 min. The solution is filtered, using a syringe filter, before injection. No additional cleanup of the sample solution is performed.
- 2. Shrimp, 1 g of shrimp is defrosted and mixed in a blender. To the 1 g of the mixed shrimp, 3 mL of water and 25 μL 10 ppb IS are added. The portion is centrifuged for 5 min (8,000 rpm). The supernatant is loaded on the cartridge and allowed to stand for 5 min. Elution is performed with 5 mL ethyl acetate. The eluate is collected and the solvent evaporated under a nitrogen stream at 40 °C. The residue is redissolved in 1 mL methanol and put in an ultrasonic bath for 1 min; the solution is filtered before injection.
- 3. Chicken, 1 g of chicken is defrosted and mixed in a blender. To the 1 g of the mixed chicken, 3 mL of water and 25 μL 10 ppb IS are added. The portion is centrifuged for 5 min (8,000 rpm). The supernatant is loaded on the cartridge and allowed to stand for 5 min. Elution is performed with 5 mL ethyl acetate. The eluate is collected and the solvent evaporated under a nitrogen stream at 40 °C. The residue is redissolved in 1 mL methanol and put in an ultrasonic bath for 1 min; the solution is filtered before injection.

#### **MRM Setting**


Name	Precursor ion	Product ion	Frag (V)	CE (V)	Dwell (ms)
TAP	354.1	185.1*	120	20	60
	354.1	289.9	120	10	60
FF	356	185.1*	120	20	60
	356	335.8	120	5	60
CAP	321.2	152.1*	120	10	60
	321.2	257.1	120	5	60
D5-CAP (ISTD)	326.2	157.2	130	15	60

#### **Results**

Name	Linearity (0.5–20 ppb)
TAP FF	0.994 0.992
CAP	0.994

#### **Sensitivity**

#### 1. CAP: 0.5 ppb S/N = 81

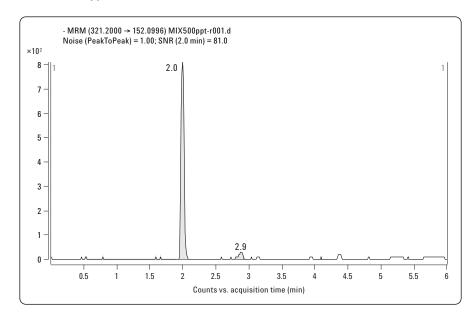


Xiaorong Ran, Tao Bo are application chemist based at Agilent Technologies, Beijing, China


### **For More Information**

For more information on our products and services, visit our Web site at www.agilent.com/chem.

#### 2. FF: 0.5 ppb S/N = 56

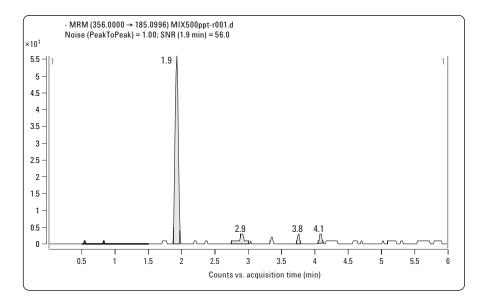

#### 3. TAP 1 ppb, S/N = 12

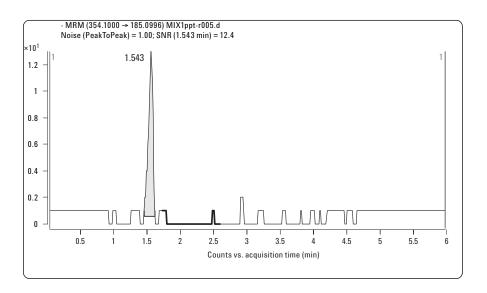

#### www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc., 2009 Printed in the USA June 15, 2009 5990-3734EN

