

# Fractionation of Formaldehyde Resin on Agilent PLgel by Gel Permeation Chromatography

# **Application Note**

Materials Testing and Research

# **Author**

Graham Cleaver
Agilent Technologies, Inc.

# Introduction

Resolving closely eluting fractions in highly disproportionate concentrations is achieved by using a heart-cutting technique. The same technique can also be used to avoid cross-fraction contamination by removing the sample matrix. In this example, heart-cutting was employed in an identification of fractions of a phenanthrene-formaldehyde resin. The resin fractions were separated by gel permeation chromatography with Agilent PLgel 10  $\mu m$  500Å columns.

# Fractionation of Phenanthrene-Formaldehyde Resin

Figure 1 shows the phenanthrene-formaldehyde resin monomer, dimer and trimer fractions. After separation of the fractions (Figure 2), identification was done using mass spectroscopy and infra-red spectroscopy.

### **Conditions**

Column Agilent PLgel 10 µm 500Å, 25 × 300 mm (p/n PL1210-6125)

Eluent Dichloromethane

Flow rate 9.0 mL/min

Conc 10%
Inj vol 2 mL
Detector RI

System Agilent PL-GPC 50





Figure 1. Structure of three fractions of a phenanthrene-formaldehyde resin.

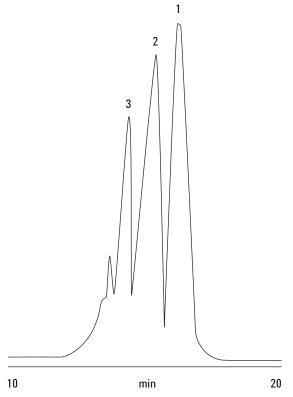


Figure 2. Separation of three fractions of phenanthrene-formaldehyde on Agilent PLgel 10 μm columns.

# **Conclusions**

Preparative gel permeation chromatography employing a heart-cut technique can be used to separate and isolate components of a polymer on the basis of size in solution.

# For More Information

These data represent typical results. For more information on our products and services, visit our Web site at www.agilent.com/chem.

# www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc., 2015 Printed in the USA April 30, 2015 5990-8603EN

